



**Catalog No. 64-030EX**

## Anti-Pertussis Toxin antibody, rabbit antiserum

Perrtussis toxin (PT) is a protein-based AB<sub>5</sub>-type exotoxin produced by *Bordeterra pertussis*. PT catalyzes the ADP-ribosylation of the  $\alpha$  subunits of the heterotrimeric guanine nucleotide regulatory proteins Gi, Go, and Gt and prevents intracellular signal transduction involving the G proteins. PT consists of one moplecule of each S1 (26 kDa), S2 (22 kDa), S3 (22 kDa), S5 (12 kDa) and two molecule of S4 (12 kDa). This product was highly purified (>90% pure) from *Bordetella pertussis* strain Tohama by the method of Skelton & Wong<sup>1</sup>. Cytotoxicity of the PT was confirmed by morphological alteration of CHO cells after treatment with 0.1 ng/ml of PT (see the Figure below).

### Applications:

1. Western blotting (1/2,000~1/10,000 dilution)
2. ELISA (1/10,000~1/20,000 dilution)
3. Dot blotting (1/2,000~1/10,000 dilution)
4. Immunoprecipitation (1/200~1/500 dilution)
5. Neutralising (Assay dependent)

Other applications have not been tested.

**Immunogen:** Immunization was Initiated with toxoid and boosted with native toxin (BioAcademia 01-503)


**Product:** Whole rabbit antiserum added with 0.09% sodium azide

**Size:** 100  $\mu$ L

**Storage:** Sent at 4°C, and upon arrival, spin-down and store at -20°C

**Data Link:** Swiss-Prot [Pertussis toxin](#)


**References:** Alouf JE & Popoff MR (Ed.) The comprehensive Sourcebook of BacterialProtein Toxins 3rd Ed. Academic Press (2006)

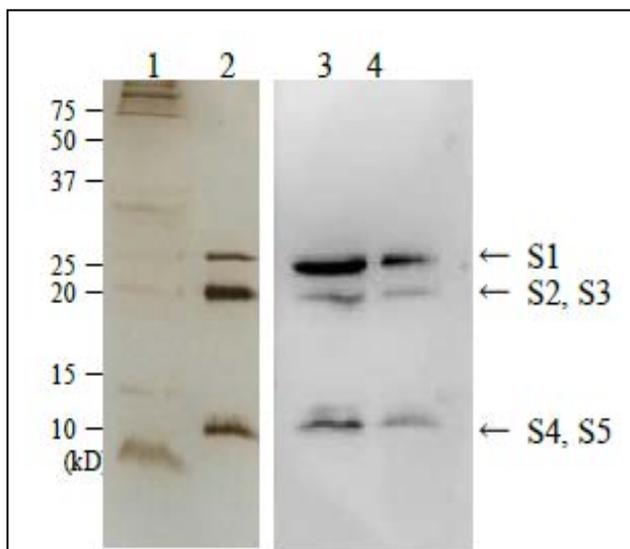

**Fig.1. Detection of perussis toxin in culture medium of Bordeterra pertussis strain Tohama by Western blotting using anti-perussis toxin antibody.**


1. Culture medium of Bordeterra pertussis. SDS-PAGE, silver-stained
2. Purified pertussis toxin (200 ng). SDS-PAGE, silver-stained
3. Western blot of culture medium of Bordeterra pertussis as in 1.
4. Western blot of purified pertussis toxin (10 ng)

The toxin consists of five subunits as indicated by S1 to S5.

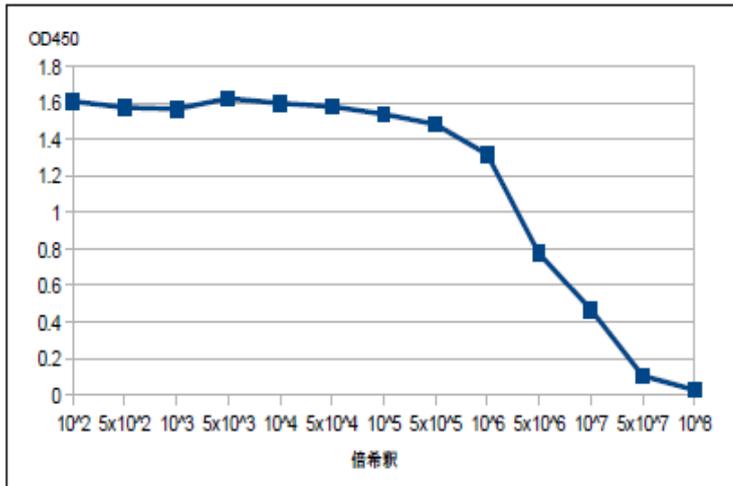

**Fig.2. Titration of antibody reactivity of anti-Pertussis antiserum by direct ELISA**

Plate was coated with 100 µg of pertussis toxin per well and 100 µl of the antiserum at the indicated dilution was added to each well and incubated. After washing, goat anti-rabbit-IgG conjugated with HRP was added as 2<sup>nd</sup> antibody. Color was developed with TMB as substrate.

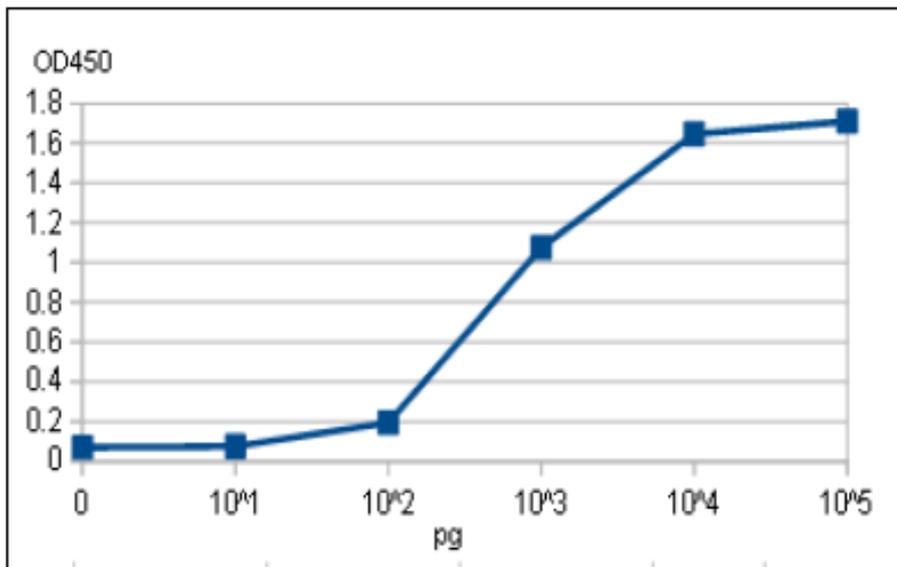

**Fig.3. Titration of pertussis toxin by direct ELISA using anti-pertussis toxin antiserum**

ELISA plate is coated with indicated amounts of pertussis toxin per well. Antiserum was used at 1/12,500 dilution. ELISA was performed as in Fig.2. Dynamic range was 100 pg to 10 ng under these conditions.

*For research use only. Not for clinical diagnosis.*

Manufactured by BioAcademia Inc.



COSMO BIO Co., LTD.  
Inspiration for Life Science

TOYO 2CHOME, KOTO-KU, TOKYO, 135-0016, JAPAN

URL: <http://www.cosmobio.co.jp>

e-mail: [export@cosmobio.co.jp](mailto:export@cosmobio.co.jp)

[Outside Japan] Phone : +81-3-5632-9617

[国内連絡先] Phone : +81-3-5632-9610

FAX : +81-3-5632-9618

FAX : +81-3-5632-9619