

Inhibitor

For research use only. Not for clinical diagnosis.

Catalog No. KN-NIGU-M02

Rat IL-1R-Ig heterodimer-containing medium for IL-1 inhibition

Background

Interleukin-I (IL-Iα and IL-1β) is a proinflammatory cytokine involved in immune responses including both innate and acquired immunity. IL-1 is thought to play a role in many diseases, including arthritis, heart disease, pancreatitis, multiple myeloma, and stroke. IL-1 receptor (IL1R) I, also known as CD121a, is an 80 kDa type I transmembrane (TM) protein that binds cytokines IL1α and IL1β and transduces a signal. Whereas IL1RII, also known as CD121b, is an 65 kDa protein that binds cytokines IL1α and IL1β but does not transduce a signal. Signal transduction requires complex formation with the IL1R accessory protein (IL1R acp), another type I TM protein. Soluble IL1R acp and soluble IL-1RII is present in normal serum and soluble form of the IL-1 receptor accessory protein (acp) increases the affinity of binding of IL-1α and IL-1β to the soluble IL1RII 1). Inhibition of IL-1 is beneficial in many animal models of disease and is expected to offer a new therapy for various human

1) Smith DE, Hanna R, Della F, Moore H, Chen H, Farese AM, MacVittie TJ, Virca GD and Sims JE: The soluble form of IL-1 receptor accessory protein enhances the ability of soluble type II IL-1 receptor to inhibit IL-1 action. Immunity 18: 87-96, 2003.

Product type Source

Recombinant Protein / Inhibitor of Rat IL-1

Rat GRO/CINC2 N (Met 1 –Asp 47)	Rat IL1R acp (Ser 21 – Tyr 356)	AAA	Rat IgG1 C region (Val 98 –Lys 326)	AQDFVQWLMNT
			S S S	
	Rat IL1R II (Met 1 – Ser 356)	AAA	Rat IgG1 C region (Val 98 –Lys 326)	Rat GRO/CINC2 C (Lys 89 – Ser 101)

Condition medium (RPMI 1640, 10%FBS, 100 unit/ml penicillin, 100µg/ml streptomycin) was obtained from Cos-7 cells cotransfected with pCAGGS- Rat GRO/CINC2αN- Rat IL-1R acp- Rat IgG1Fc and pCAGGS- Rat IL-1R II- Rat IgG1 Fc- Rat GRO/CINC2αC by Fugene 6 (Roche, Indianapolis, IN). It was filtered by Millex-GV Filter Unit, 0.22 μm (Millipore, Billerica, MA). It contains not only Rat IL-1R-Ig heterodimer but also IL-1R acp-IgG1Fc homodimer and IL-1R II- IgG1Fc homodimer. However, IL-1 inhibition effect by IL-1R-lg heterodimer is much stronger than that by their homodimers. MOCK medium was obtained from them trasfected with pCAGGS.

Form Liquid (control medium is attachd)

Volume/ Concentration 500 μl / tube, Rat IL-1R-lg heterodimer 100 fmol/ml


Concentration of Rat IL-1R-Ig heterodimer in condition medium, which is measured by Rat

GRO/CINC2α ELISA kit (IBL, Takasaki, Japan), is approximately 100 fmol/ml.

Activity

Measured by its ability to inhibit MCP-1 gene expression in NRK-49F cell (rat kidney fibroblast) after culture for 24 hr with rrlL-1α and the condition medium. approximately 0.625 fmol/ml and 1.25 fmol/ml IL-1R-la heterodimer-containing medium will inhibit 88% and 96% of the biological response due to 200 pg/mL of rrIL-1a. ID50; 0.15 fmol/ml. MOCK medium will not inhibit it.



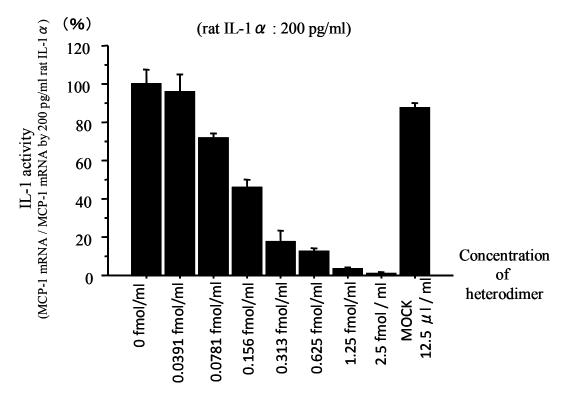

Storage

Store below -20°C (below -70°C for prolonged storage).

Aliquot to avoid cycles of freeze/thaw.

Reference

 IL-1 Receptor Accessory Protein-lg/IL-1 Receptor Type II-lg Heterodimer Inhibits IL-1 Response More Strongly than Other IL-1 Blocking Biopharmaceutical Agents. J Clin Immunol. 2010 Dec 22. (PMID: <u>21181432</u>)

For research use only. Not for clinical diagnosis.



COSMO BIO CO., LTD.

Inspiration for Life Science

TOYO 2CHOME, KOTO-KU, TOKYO, 135-0016, JAPAN

URL: http://www.cosmobio.co.jp e-mail: <u>export@cosmobio.co.jp</u> [Outside Japan] Phone: +81-3-5632-9617 [国内連絡先] Phone: +81-3-5632-9610

FAX: +81-3-5632-9618 FAX: +81-3-5632-9619