

Coat Quick “G”

Catalog No. DAI-CQG

Principle

Coat Quick "G" Pen provides a thin layer of adherent coating on the grid. The grid will be ready for tissue section mounting after drying for one or two minutes at room temperature.

Coat Quick "G" Pen is also effective as a pretreatment prior to application of supporting films such as former, collodion and other delicate membranes by minimizing any dislodging or widening of a break in the supporting film. This procedure is very easy and time saving, since the coated grid will not stick to the tip of the pen or to filter paper. This is an excellent adjunct for specimen retrieval techniques with TEM grid.

How to Use

1. Press the tip of the Coat-Quick "G" Pen straight down several times or until the fluid starts to soak the tip.
2. A light touch with this pen will provide a thin layer of adhesive on the grid. Layer will dry in 1-2 minutes at room temperature and is then ready to receive the sections.
3. The grid will be ready for tissue section mounting.
4. Keep tightly capped after using.

It contains 2 pieces.

Store in a room temperature



COSMO BIO CO., LTD.
Inspiration for Life Science

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME, KOTO-KU, TOKYO 135-0016 JAPAN
TEL: (81)3-5632-9617 / FAX: (81)3-5632-9618 / e-mail: export@cosmobio.co.jp / URL: www.cosmobio.com