PAP Pen Super-Liquid Blocker

This ready-to-use hydrophobic barrier pen is designed for immunohistochemistry applications. The surface tension provided by the circle, square, or any other shapes drawn to fit tissue sections with the pen, ensures that only the amount of antibody solution needed for the sufficient reaction will be applied. The pen is stable up to 129 degrees C in microwave heat.

The PAP Pen is an effective tool for immunostaining procedures by the Peroxidase-Antiperoxidase (PAP) method, ABC method, LSAB method, polymeric method, immunofluorescence method, ASD method, Enzyme method and Frozen Section method.

For paraffin sections, the PAP Pen should be applied after deparaffinization. For frozen sections, the PAP Pen can be applied after fixation.

Directions to Use:

- Shake well before use. Push down on the tip a couple times to release liquid into the tip.
- Deparaffinize tissue sections and hydrate to water. Perform antigen retrieval if necessary.
- Wipe away excess liquid around the section on slide with Kimwipes or tissue paper.
- Allow tip to get moderately wet. Encircle the tissue sections or draw lines on both sides of the section and let dry for 10 to 15 seconds.
- Apply drops of washing buffer to cover the section.
- Proceed with standard immunohistochemistry procedure.

Size and Specification

- Content: 6cc
- Size: 141mm(L) x 15mm(D)
- Pen Tip: 4mm

For research use only, not for diagnostic use

COSMO BIO Co., LTD.

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU. TOKYO 135-0016, JAPAN
TEL : (81)3-5632-9617
FAX : (81)3-5632-9618
e-mail : export@cosmobio.co.jp
URL : www.cosmobio.com


Precautions

The push-and-flow method has to be applied for stationary with fluids that may separate,. If the pressure inside the pen rises, it may lead to leakage of the fluid.

Before using the pen, please make sure you do what is shown in the figure "How to release air." Compared with regular stationary equipment, the PAP pen uses a chemical that vaporizes easily. When using the pen after long periods of storage or when using it for the first time, please carry out the "air-release" procedures. Also, please note that if too much pressure is applied when using the pen, the closing valve will open and more fluid will come out.

Ways to avoid fluid leakage of PAP-Pen

1. Before use, direct the tip of the Pen to the ceiling and push the tip down and release air. #1
2. Direct the tip of the Pen down, and wait until the fluid comes out of the pen. #2 If the pen is shaken up and down, fluid will accumulate in the tip holder and can cause fluid leakage.
3. Please carry out #1 again in order to allow the excess fluid in the tip holder to return to the body of the Pen.
4. If no fluid comes out, please repeat #1 to #3.

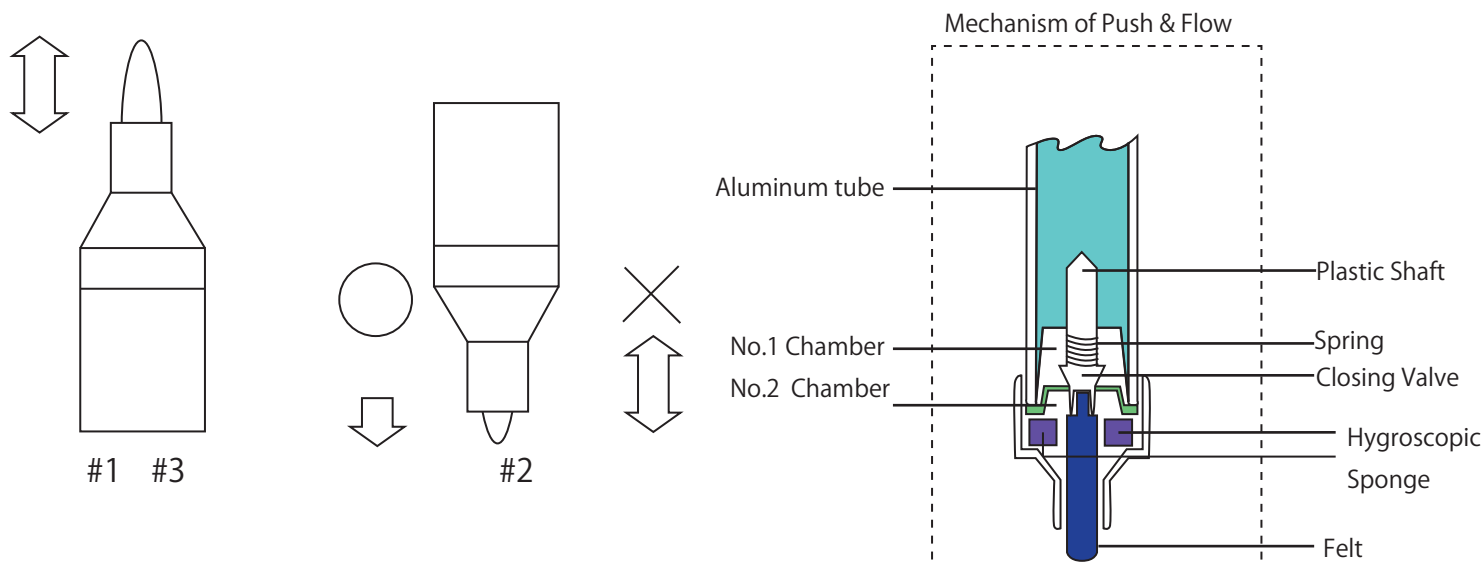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COSMO BIO CO., LTD.
DAIDO SANGYO CO., LTD.