



Anti Estrone-3-Glucuronide polyclonal antibody

Product No. FKA-224

The Cross reactivities of anti Estrone-3-Glucuronide (FKA-224) by RI Assay

Antigen	Estrone-3-Glucuronide-BSA
Antibody[FKA-224]	lyophilized Serum 2.5μl / Vial
Dilution of Serum	2000test for RI Assay 250μl / tube use
Label	Estrone[2,4,6,7-3H] 5000dpm/tube use
Binding of Total	about 50%
Estrone dose of B0/2	200 pg
Strage/Handling	Maintain at 4°C-8 °C (-20 °C for long term storage) in undiluted aliquots. Avoid repeated freeze/thaw cycles.

Steroid	Cross Reaction
Estrone (E1)	100 %
Estrone(E1)-3-Sulfate	40.00 %
Estrone(E1)-3-Glucuronide	170.60 %
Estradiol(E2)	1.00 %
Estradiol(E2)-3-Sulfate	5.00 %
Estradiol(E2)-3-Glucuronide	1.50 %
Estradiol(E2)-3-Sulfate	0.00 %
Estradiol(E2)-17-Glucuronide	0.10 %
Estriol(E3)	0.00 %
Estriol(E3)-16-Glucuronide	0.00 %
Estriol(E3)-Sulfate	0.00 %
Testosterone	0.00 %
The other steroids	0.00 %

For research use only; not for use as a diagnostic.

