



Code No. KAL-KH036-EX

For research use only

Anti Nitrotyrosin Monoclonal Antibody

(Clone No. 2H1)

To elucidate the function of Nitric Oxide (NO) related signal transduction, we developed new monoclonal antibody to Nitrotyrosin (Clone No.2H1). There are two pathways, which is engaged in the signal transduction regarding vascular relaxation on endothelial cell. One is through activated guanylate cyclase that is cGMP dependent and another is cGMP independent pathway which requires reactive NO derived substance such as, peroxynitrate(ONOO^-), N_2O_3 , $\text{N}_2\text{O}_4(\text{NO}_2)$.

Recently, various functions of reactive NO derived substance has been identified for example regulating protein phosphorylation and inducing apoptosis. So, it is believed that reactive NO derived substance is involved in signal transduction among cells. Immunohistochemical study revealed that nitotyrosin residue is produced in some patient such as athrosclerosis, Alzheimer's disease, Parkinson's disease, and acute lung damage.

This antibody is very useful for the research of reactive NO derived substance.

Package Size	100 μ g (400 μ L / vial)
Format	Mouse monoclonal antibody 0.25mg/mL
Buffer	Block Ace as a stabilizer, containing 0.1% Proclin as a bacteriostat
Storage	Store below -20°C Once thawed, store at 4°C . Repeated freeze-thaw cycles should be avoided
Clone No.	2H1
Subclass	IgG1
Purification method	The spleen cells from BALB/c mouse, immunized with nitrotyrosin-HSA, were fused to myeloma P3U1 cells. The screening of the hybridoma cells was performed on ELISA. The cell line was grown on non-serum medium, from which the antibody was purified by Protein G affinity chromatography.



For research use only

Anti Nitrotyrosin Monoclonal Antibody

(Clone No. 2H1)

【Reference】

1. Furchgott, R. F. *Acta Physiol. Scand* 139, 257-70 ,1990
2. Moncada, S. &Higgs, A. *N. Eng. J. Med* 329, 2002-12 ,1993
3. Ischiropoulos, H. *Arch. Biochem. Biophys* 356, 1-11 ,1998
4. Stamler, J. S., Toone, E. J., Lipton, S. A. & Sucher, N. J. *Neuron* 18, 691-6 ,1997
5. Akaike, T., et al. *J. biochem* 122, 459-66 ,1997
6. Beckman, J. S., Ye, Y. Z., et al. : Extensive nitration of protein tyrosines in human atherosclerosis detection by immunohistochemistry. *Biol. Chem. Hoppe-Seyler* 375, 81-88, 1994
7. J. S. Luoma, P. Stralin, et al. : Expression of extracellular SOD and iNOS in macrophages and smooth muscle cells in human and rabbit atherosclerotic lesions. Colocalization with epitopes characteristic of oxidized LDL and peroxynitrite-modified proteins. *Atheroscler. Thromb. Vasc. Biol* 18, 157-167, 1998
8. P. F. Good, P. Werner, et al. : Evidence for neuronal oxidative damage in Alzheimer's disease. *Am. J. Pathol* 149, 21-28, 1996
9. P. F. Good, A. Hsu, et al. : Protein nitration in Parkinson's disease. *Journal of Neurophthology and Experimental Neurology* 57, 338-342, 1997
10. I.G.Haddad, G. Pataki, et al. :Quantitation of nitrotyrosine levels in lung section of patients and animals with acute lung injury. *J. Clin. Invest* 94,2407-2413,1994

* These references are the background of nitrotyrosin , and are not this antibody examples .

Distributor



COSMO BIO CO., LTD.
Inspiration for Life Science

TOYO EKIMAE BLDG. 2-20, TOYO 2CHOME
KOTO-KU, TOKYO 135-0016, JAPAN
TEL : +81-3-5632-9617
FAX : +81-3-5632-9618
URL : <http://www.cosmobio.co.jp/>
e-mail : export@cosmobio.co.jp

Manufacturer

Trans Genic Inc.

7-1-6 Minatojimmaminami-machi,
Chuo-ku, Kobe,650-0047 JAPAN
TEL : +81-78-306-0590
FAX : +81-78-306-0589
URL : <http://www.transgenic.co.jp/>
e-mail : techstaff@transgenic.co.jp