

# Cryopreservation of Mouse Spermatozoa in *FERTIUP*<sup>®</sup>

Department of Reproductive Engineering, Center for Animal Resources and Development, Kumamoto University, Japan  
Senior Editor: Naomi Nakagata

Center for  
Animal  
Resources and  
Development  
20140415


## -- Materials and Equipment --

1. Male mice (12-70 weeks old)
2. Micro-spring scissors (5 mm blade)
3. Pair of watchmaker's #5 forceps
4. FERTIUP<sup>®</sup> (Cryoprotectant : CPA ; COSMOBIO Co., LTD. , Cat. no. KYD-001-EX)
5. HTF (COSMOBIO Co., LTD. , Cat. no. CSR-R-B070)
6. Plastic dish (35mm X 10mm Cat.No.430588; CORNING)
7. Pipette tips
8. 0.25mL PETG cotton-plugged Sperm Straws (COSMOBIO Co., LTD. , Cat. no. KYD-S020X10)
9. Micropipettes
10. Straw connector (COSMOBIO Co., LTD. , Cat. no. KYD-S025)
11. Impulse sealer
12. Freezing canister (COSMOBIO Co., LTD. , Cat. no. KYD-S018)
13. Cryobiological container
14. Hot plate

## -- Procedure --


### ●●Preparation the Freezing Canister●●

- 1 Insert a piece of styrofoam tightly into the bottom of the syringe.
- 2 Heat seal the mouth of the syringe tip.
- 3 Fix the syringe to the acrylic bar.

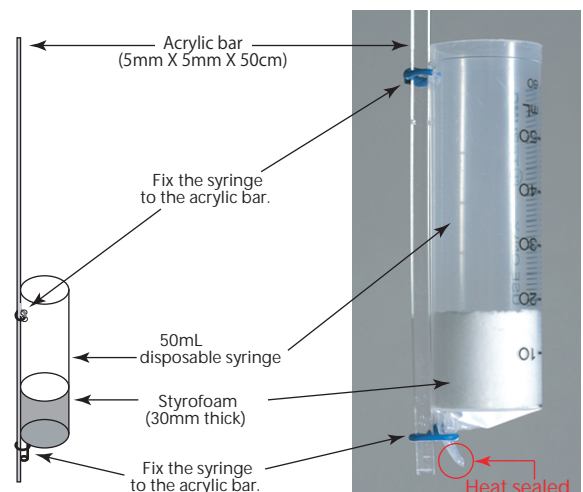


### ●●Preparing a Straw Connector●●

- 1 Using a 1mL syringe, a 3-way stopcock, a piece of silicone tube, a plastic tube and a silicone cap, make a straw connector.
- 2 To use the straw connector, cut away the cotton plug from a straw (0.25mL French Straw), then attach the straw to the silicone cap at the end of the connector.

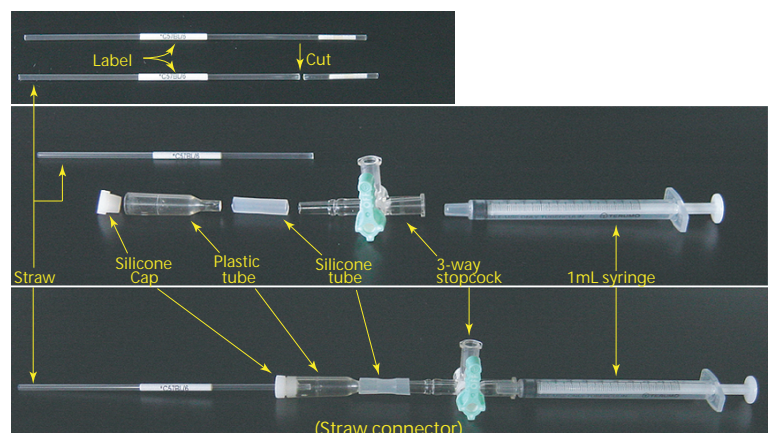


## ●●Preparing Sperm Suspension●●

- 1 Make a drop of 60  $\mu\text{L}$  of FERTIUP<sup>®</sup> (CPA) on a 35mm culture dish and cover with liquid paraffin.
- 2 Add a 60  $\mu\text{L}$  aliquot of the same solution to the drop (final volume: 120  $\mu\text{L}$ ) to make a tall, semispherical drop.
- 3 Sacrifice a male mouse (>12 weeks-old) via cervical dislocation and remove the two cauda epididymides aseptically.
- 4 Place the cauda epididymides on a piece of filter paper and completely remove any fat and blood under a microscope.
- 5 Transfer the cauda epididymides into the drop of FERTIUP<sup>®</sup> (CPA) and use a pair of watchmaker's #5 forceps and micro-spring scissors to make 5 or 6 incisions in the epididymides.
- 6 Place the dish on a hot plate at 37°C for 3 minutes. During this time, rotate the dish every minute to disperse sperm from the organs in the FERTIUP<sup>®</sup> (CPA).

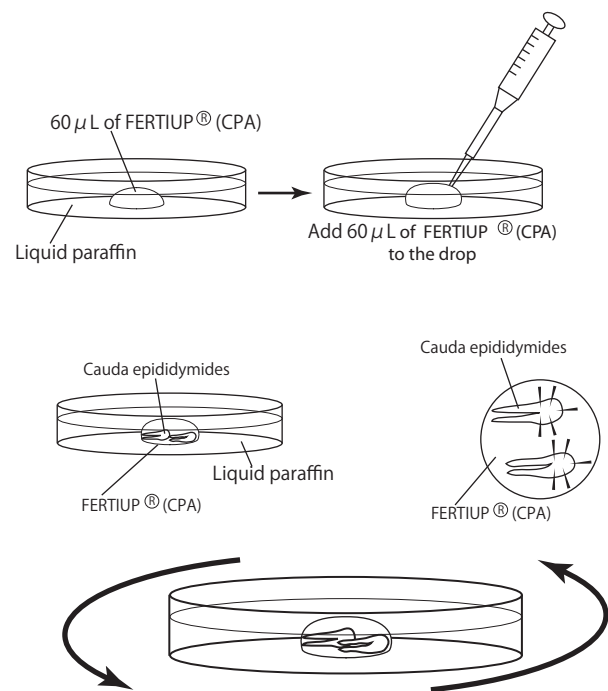

## ●●Preparing Freezing Straw Containing Sperm Suspension●●


- 1 Connect a straw to a straw connector.
- 2 Carefully aspirate the contents into the straw in following order:
  - a. 100  $\mu\text{L}$  of HTF,
  - b. 15mm of air,
  - c. 10  $\mu\text{L}$  of the sperm suspension,
  - d. Another 15 mm of air.
- 3 Seal both sides of the straw using an impulse sealer.  
 Comment: Loading 100  $\mu\text{L}$  of HTF into the straw prevents the straw from floating on the surface of liquid nitrogen. This is because the HTF acts as a weight that forces the straw to sink into the liquid nitrogen.

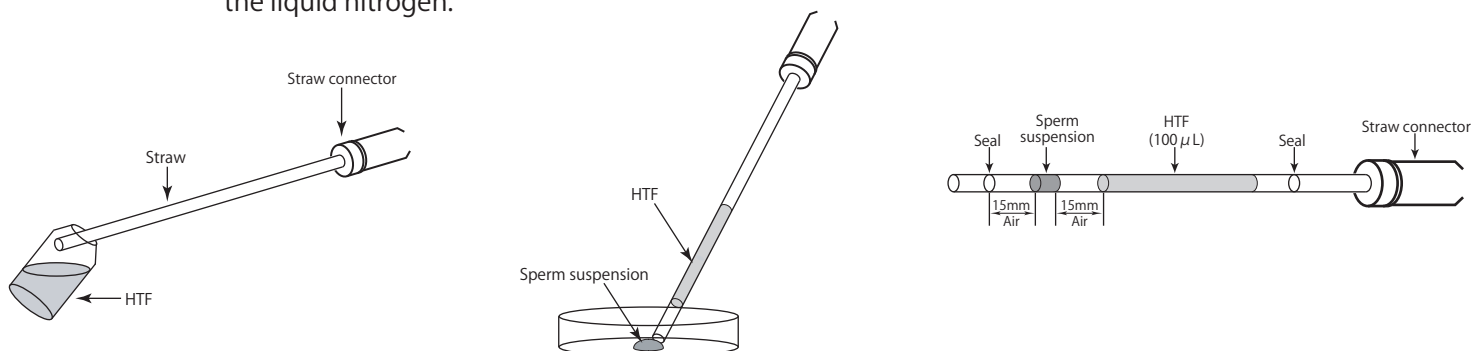

- 4 Create 10 samples per mouse in the same manner as described above.

## ●●Sperm Freezing using a Cryobiological Container●●

- 1 Put the samples into a freezing canister and float them on liquid nitrogen in a cryobiological container.
- 2 After 10 minutes, quickly immerse the freezing canister in the liquid nitrogen.

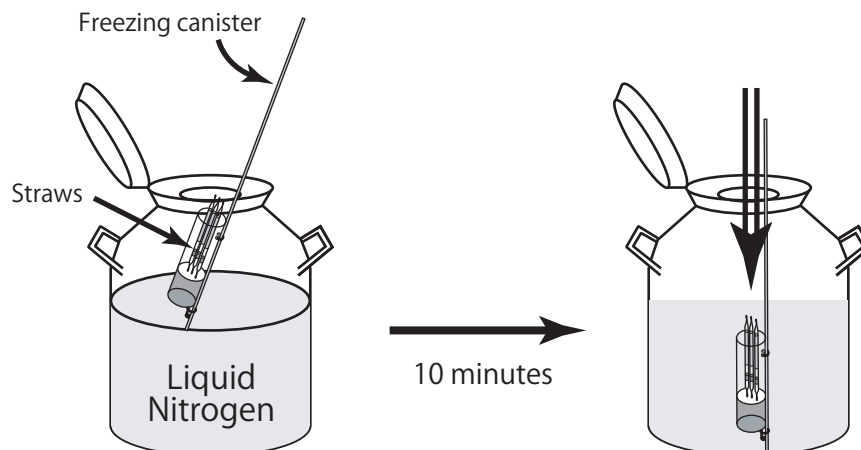

# *In Vitro* Fertilization using Spermatozoa Cryopreserved in *FERTIUP*<sup>®</sup>

Department of Reproductive Engineering, Center for Animal Resources and Development, Kumamoto University, Japan  
Senior Editor: Naomi Nakagata




## --Materials and Equipment--

1. Female mice superovulated with PMSG and hCG
2. FERTIUP<sup>®</sup> (Preincubation Medium : PM ; COSMOBIO CO.,LTD. Cat. no. KYD-002-EX )
3. CARD MEDIUM<sup>®</sup>(COSMOBIO CO.,LTD. Cat. no. KYD-003-EX)
4. HTF (COSMOBIO Co., LTD. , Cat. no. CSR-R-B070)
5. Pipette tips (Pipette Tip Cat.No.114 Quality Scientific Plastics)
6. Plastic dishes ( 35mm X 10mm Cat.No.430588; CORNING)
7. Straw connector (COSMOBIO CO.,LTD. Cat. no. KYD-S025)
8. Water bath maintained at 37°C
9. Float for thawing
10. Micropipettes
11. Humidified incubator (37°C, 5% CO<sub>2</sub>, 95% air)

## --Procedure--

### ●●Preparation of the Float for Thawing●●

- 1 Using some styrofoam and a 50mL plastic centrifuge tube, make the float as shown in the diagram below.

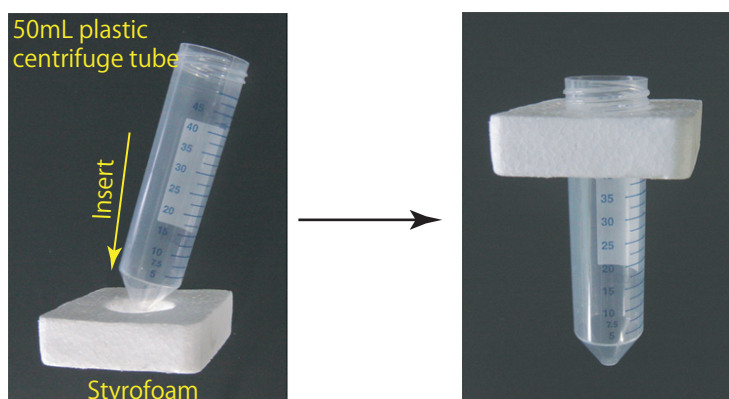


### ●●Preparation for Thawing●●

- 1 Prepare a water bath maintained at 37°C.
- 2 Pour water (37°C) into the 50mL plastic centrifuge tube part of the styrofoam/centrifuge tube assembly, and float it in a water bath.
- 3 Put 1 drop (90  $\mu$ L / drop) of FERTIUP (PM)<sup>®</sup> into a dish and cover it with liquid paraffin 30 minutes before thawing a frozen straw, and place the dish in an incubator (37°C, 5% CO<sub>2</sub> in air).


## ●●Thawing the Mouse Spermatozoa●●

- 1 Remove a frozen straw from the liquid nitrogen and hold it in the air for 5 seconds.
  - 2 After completing step 1, immediately immerse the frozen straw in the styrofoam/centrifuge tube assembly (in a water bath maintained at 37°C) for 10 minutes.
- Note:** To ensure warming of the frozen sperm, completely immerse the part of the straw containing the sperm in the water bath. Furthermore, frozen-thawed mouse spermatozoa are sensitive to environmental changes. If the straw is not kept in the water bath long enough (10 minutes), the motility of the cryopreserved spermatozoa will be reduced.
- 3 10 minutes after immersion, remove the straw from the styrofoam/centrifuge tube assembly.
  - 4 Use fine tissues to wipe any water from the straw.

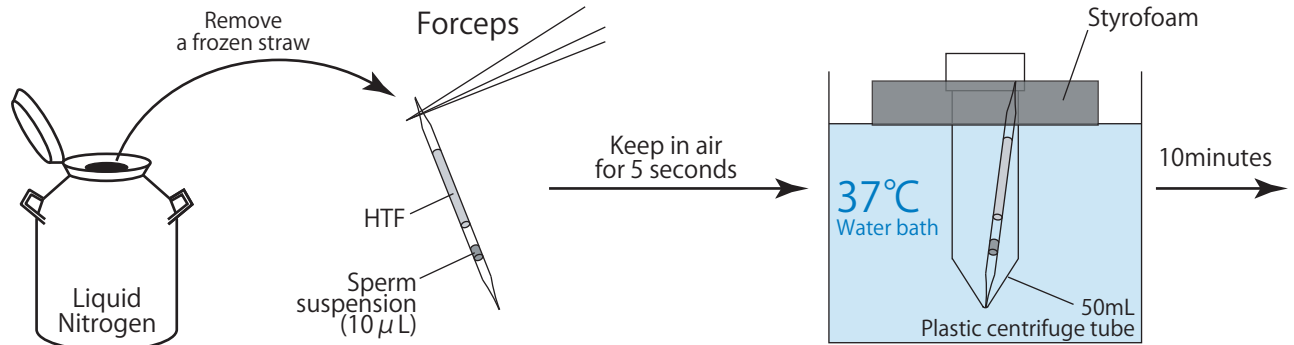


## ●●Transferring and Preincubating the Thawed Sperm Suspension●●

- 1 Cut the sealed end of the straw furthest from the sperm and attach the cut end to the straw connector (see figure below). Then cut the other sealed end from the straw.

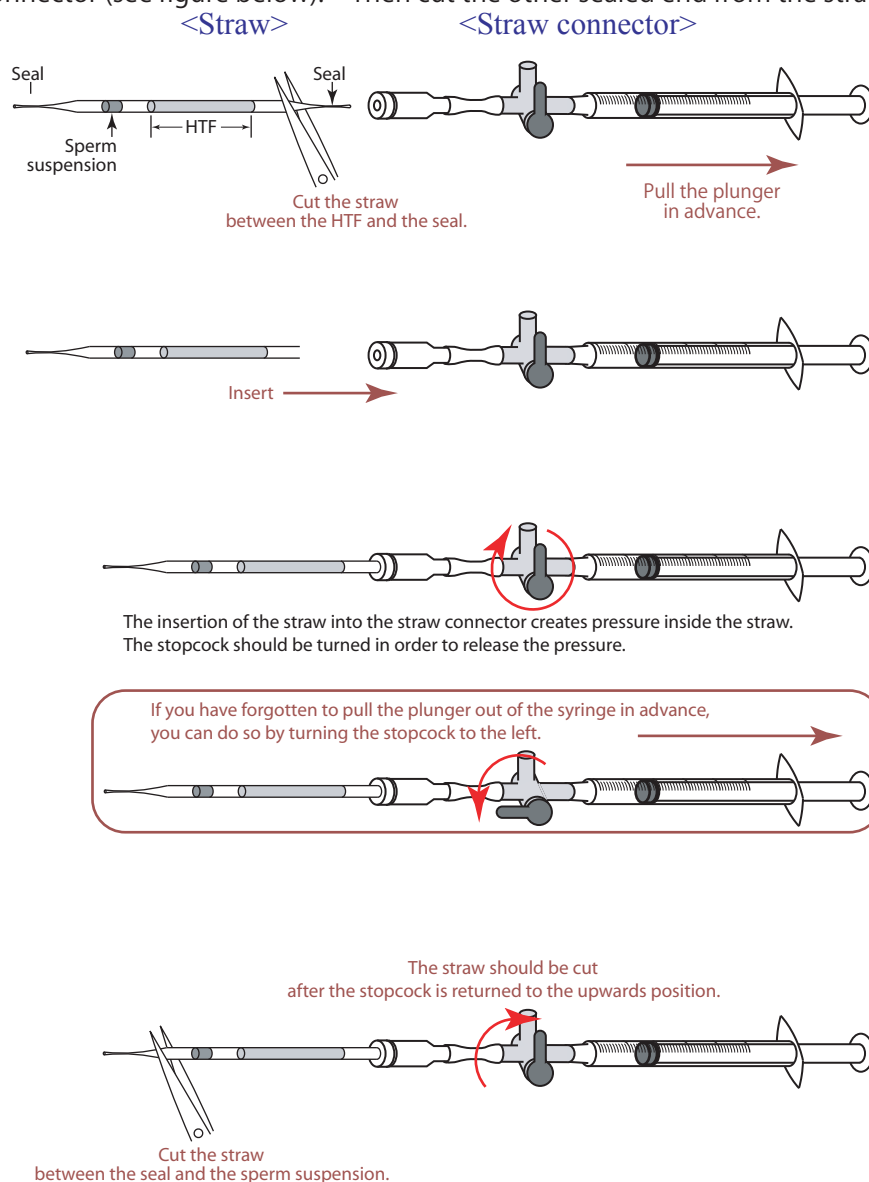

- 2 Push the plunger to transfer only the sperm suspension into the drop of FERTIUP<sup>®</sup> (PM), and place the dish in an incubator (37°C, 5% CO<sub>2</sub> in air) for 30 minutes.

**Note:** Do not disturb the dishes containing cryopreserved spermatozoa until they are moving sufficiently within the medium. If the dishes are disturbed before the spermatozoa start to move, then they will not recover full motility.

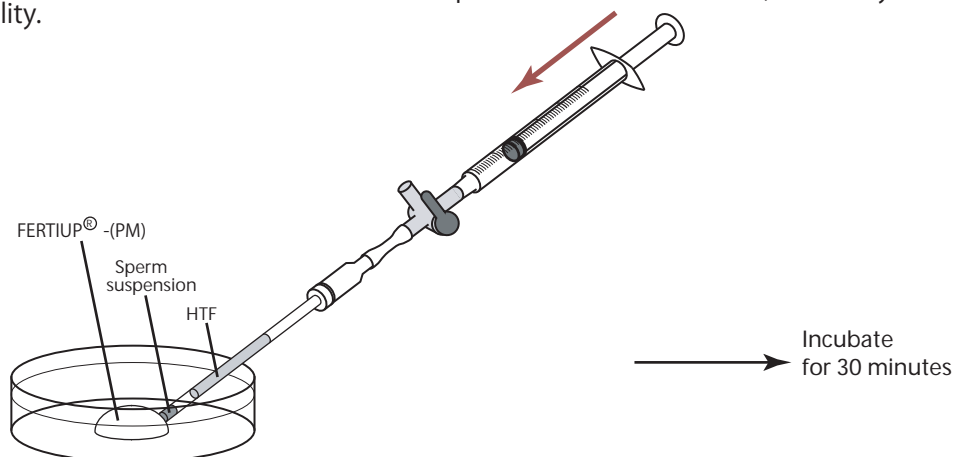

## ●●Preparation of Dishes for Fertilization and Washing●●

- 1 Put 1 drop (90  $\mu$ L / drop) of CARD MEDIUM<sup>®</sup> into a dish and cover it with liquid paraffin 10 mins before collecting of oocytes, and place the dish in an incubator (37°C, 5% CO<sub>2</sub> in air).
- 2 Put 4 drops (80  $\mu$ L / drop) of HTF into a dish and cover them with liquid paraffin. Place the dish in an incubator (37°C, 5% CO<sub>2</sub> in air) for at least 30 minutes.

## ●●Insemination●●

- 1 Sacrifice female mice 15-17 hours after an hCG injection and remove the oviducts.
- 2 Using fine, sharp needles, release up to 4-6 cumulus-oocytes-complexes (COCs) masses into each drop of CARD MEDIUM<sup>®</sup> (90  $\mu$ L).

**Note:** Be sure to carry out all operations, from sacrificing the female and removing her oviducts to introducing the COCs into a drop of CARD MEDIUM<sup>®</sup>, in the shortest time possible (within 30 seconds). Moreover, when carrying out this process alone, do not sacrifice multiple mice at once; instead, sacrifice one mouse and swiftly remove its oviducts before moving on to the next mouse.


- 3 Using a wedge-shaped pipette tip (Pipette Tip Cat.No.114 Quality Scientific Plastics), aspirate 10  $\mu$ L of the preincubated sperm suspension from the edge of the drop.

**Comment:** Spermatozoa with high motility have a tendency to gather near the edge of the drop.  
**Comment:** It is possible to aspirate 10  $\mu$ L of sperm suspension 3-4 times per drop.


- 4 Add 10  $\mu$ L of sperm to each drop of fertilizing CARD MEDIUM<sup>®</sup> containing the COCs.

**Note:** Perform the pipette operation mentioned in steps 3 and 4 as gently as possible.

- 5 Incubate the oocytes and spermatozoa for 3 hours in an incubator (37°C, 5% CO<sub>2</sub> in air).

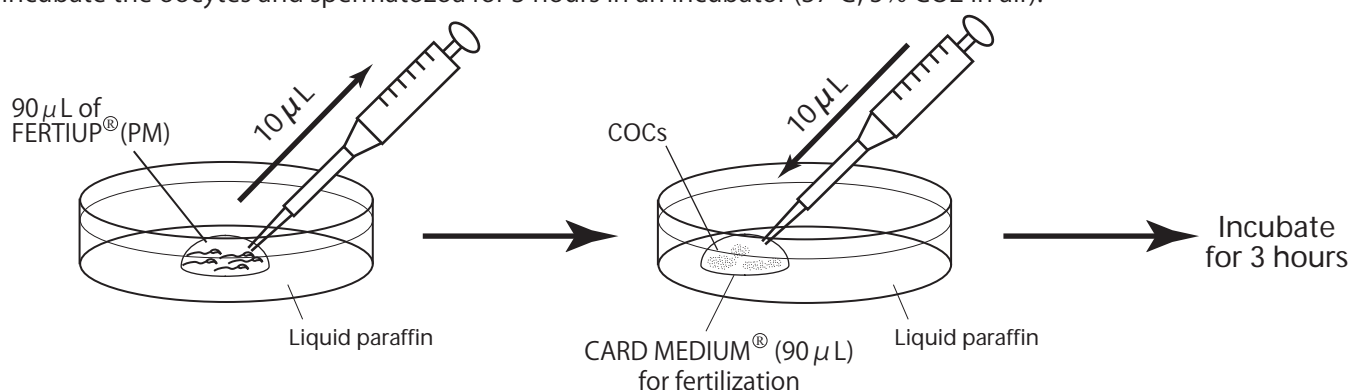


[Aspirating the Sperm Suspension from the Edge of the Drop]

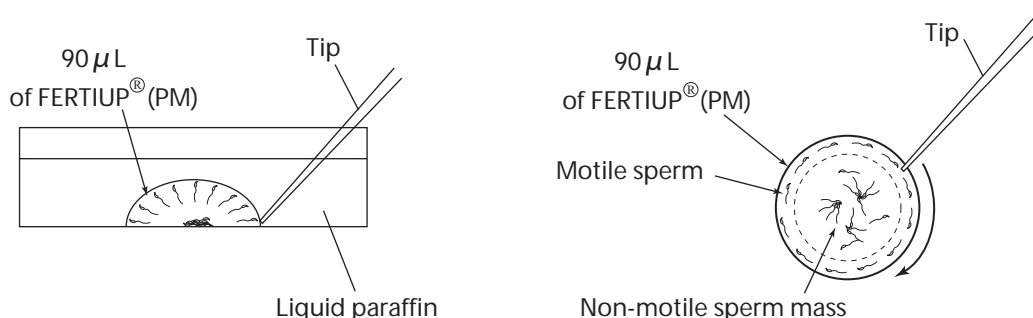




- 6 After incubating for 3 hours, wash the oocytes 3 times in fresh HTF (80  $\mu$ L) in a washing dish, avoiding transfer of CARD MEDIUM®.

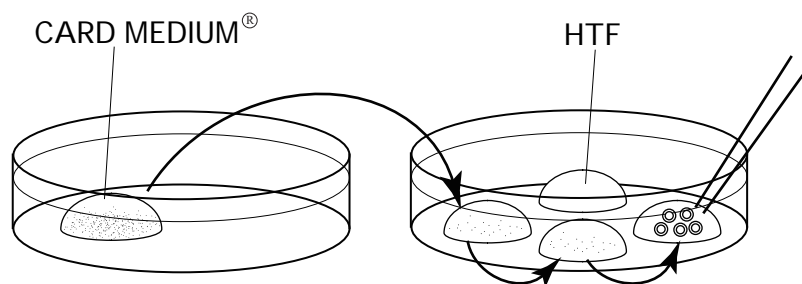

- 7 6 hours after insemination, observe them in the third drop of HTF and remove any parthenogenetic oocytes which have only one pronucleus.
- 8 After overnight culture of the oocytes, transfer the obtained 2-cell stage embryos only to the fourth drop of HTF. These embryos can now be vitrified or transferred.

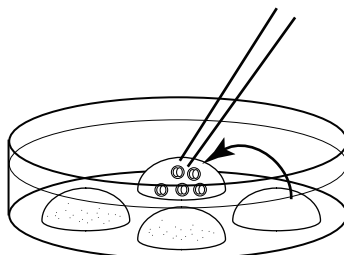

## —References—

For information on cryopreservation techniques for mouse spermatozoa, or on techniques used in thawing, *in vitro* fertilization and obtaining pups through embryo transfer, please refer to CARD website below.  
<http://card.medic.kumamoto-u.ac.jp/card/english/sign/index.html>

### ●●References●●

- 1) Takeo T., Hoshii T., Kondo Y., Toyodome H., Arima H., Yamamura KI., Irie T., and Nakagata N. 2008. Methyl-beta-cyclodextrin improves fertilizing ability of C57BL/6 mouse sperm after freezing and thawing by facilitating cholesterol efflux from the cells. *Biol Reprod.* 78(3): 546-51.
- 2) Takeo T., and Nakagata N. 2010. Combination medium of cryoprotective agents containing L-glutamine and methyl-  $\beta$ -cyclodextrin in a preincubation medium yields a high fertilization rate for cryopreserved C57BL/6J mouse sperm. *Lab Anim.* 44(2): 132-7.

## —Note—

For research use only. Not for human or medicinal use.

FERTIUP Cryoprotectant may precipitate when it is stored at low temperature.  
 In case the precipitant is obtained, please incubate the vial 30 min to 1 hour at 37°C. After dissolving, there is no problem in use.

\*FERTIUP® was jointly developed by KYUDO CO., LTD. and Kumamoto University.

\*FERTIUP® is a trademark of KYUDO CO., LTD.

\*FERTIUP® PATENT Number JP4683408

\*CARD MEDIUM® are based on the results of research in The Center for Animal Resources and Development (CARD), Kumamoto University, Japan.

\*CARD MEDIUM® is a trademark of KYUDO CO., LTD.

Manufacturer



COSMO BIO Co., LTD.

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,  
 KOTO-KU. TOKYO 135-0016, JAPAN  
 TEL : (81)3-5632-9617  
 FAX : (81)3-5632-9618  
 e-mail : export@cosmobio.co.jp  
 URL : www.cosmobio.com

## Instruction manual

CARD MEDIUM<sup>®</sup> is a novel medium which helps attain a high fertilization rate *in vitro* in mice, especially when used in combination with FERTIUP<sup>®</sup> pre-incubation medium.

CARD MEDIUM<sup>®</sup> was developed by

The Center for Animal Resources and Development (CARD), Kumamoto University, Japan.

### Contents

An ampoule including medium (A)

A 1.5mL plastic tube (C)

A 2.5mL disposable syringe

A filter unit (pore size: 0.22  $\mu$ m)

A vial including powder (B)

A 1.5mL plastic tube (D)

A needle

### Storage


Store at 4 °C. Please see the labels on the ampoule (A) and the vial (B) for the use-by date.

### Note


For research use only. Not for human or medicinal use.

### Preparation of CARD MEDIUM<sup>®</sup> for *in vitro* fertilization


- 1) Take 1mL of the medium from the ampoule (A) and add it into the vial (B). Close the lid, then mix the medium and the powder in the vial.

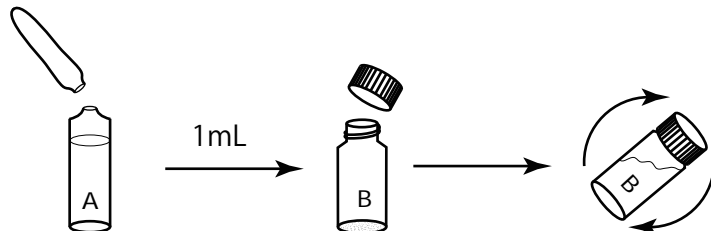

- 2) Take 1.5mL of the medium from the ampoule (A) and add it into the tube (C).

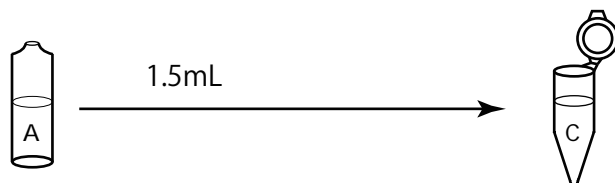


- 3) Put predetermined volume\* from the vial (B) into the tube (C) and mix them.

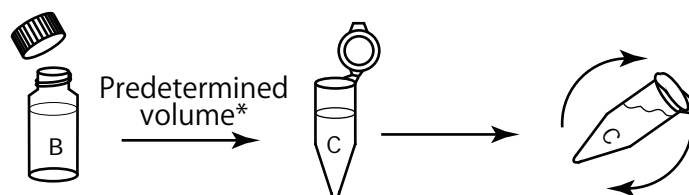


Sperm	Predetermined Volume*
Frozen-thawed	15 $\mu$ L
Cold stored	7.5 $\mu$ L
Fresh	3.75 $\mu$ L


4) Connect the needle to the syringe and suck up all of the medium in the tube (C).

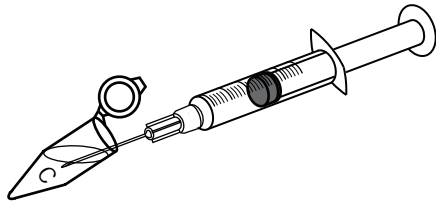

5) Remove the needle from the syringe and connect the filter unit to the syringe.

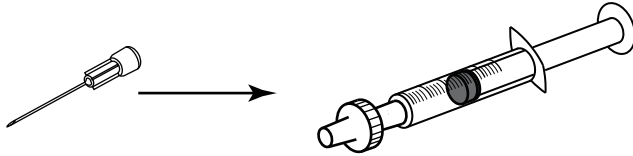


6) Pushing the syringe in, discard 2~3 drops of medium and inject the remaining medium into the tube (D).

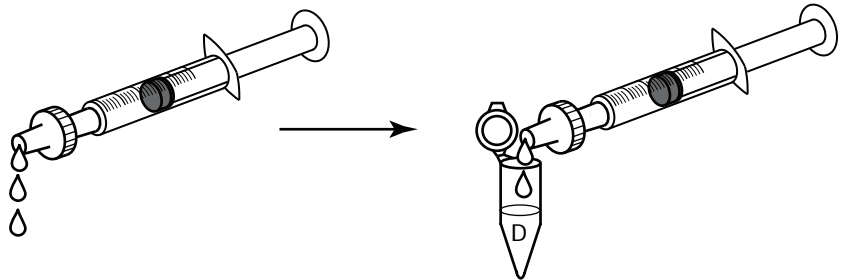

7) Make a drop\* of the medium in the tube (D) in a fertilization dish, then keep the dish in an incubator (37°C, 5% CO<sub>2</sub> in air) for 10 minutes.

\*A drop volume of CARD MEDIUM<sup>®</sup> for fresh and cold stored sperm: 200  $\mu$  L  
A drop volume of CARD MEDIUM<sup>®</sup> for frozen-thawed sperm: 90  $\mu$  L

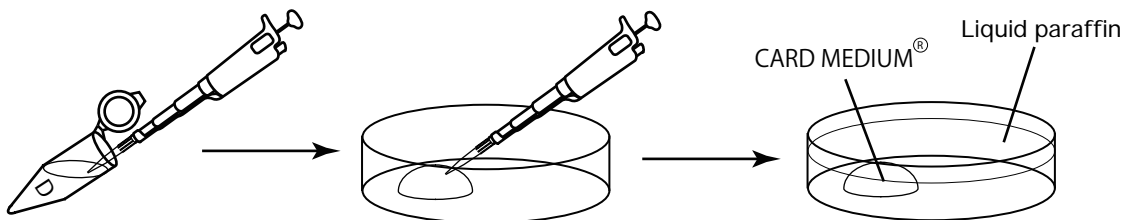

8) Introduce oocytes from superovulated female mice into the drop of CARD MEDIUM<sup>®</sup>, then add preincubated sperm into the drop to fertilize them.

**Note:** Once the ampoule (A) is opened, the medium within cannot be stored. Moreover, once the medium has been used to dissolve the powder in the vial (B), the medium cannot be stored, even if diluted or filtered. Please use the medium within 4-5 hours of step 1).

**Comment:** FERTIUP<sup>®</sup> is a recommended medium for the preincubation of sperm.

\*CARD MEDIUM<sup>®</sup> is a trademark of Kumamoto University

\*CARD MEDIUM<sup>®</sup> are based on the results of research in The Center for Animal Resources and Development (CARD), Kumamoto University, Japan.

\*FERTIUP<sup>®</sup> is a trademark of KYUDO CO., LTD.

\*FERTIUP<sup>®</sup> was jointly developed by KYUDO CO., LTD. and Kumamoto University.

\*FERTIUP<sup>®</sup> PATENT Number JP4683408

Manufacturer



COSMO BIO Co., LTD.

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,  
KOTO-KU. TOKYO 135-0016, JAPAN  
TEL : (81)3-5632-9617  
FAX : (81)3-5632-9618  
e-mail : export@cosmobio.co.jp  
URL : www.cosmobio.com