

# ALP Stabilizer—ST3010, ST4010

—Synthetic polymer type alkaline phosphatase stabilizer—

## —Features—

- ALP stabilizer for enzyme Immunoassay
- Do not contain any bio-derived materials
- Excellent stability for ALP conjugate protein




## —Usage—

- Use it like usual stabilizer
- Use from 1 to 25-fold dilution --- recommend stock solution ---
- Recommend dilute ALP conjugate over 10 fold

## —Precautions for use—

The product contain 0.09% sodium azide as a preservatives. Be careful of intake by mistake or waste.

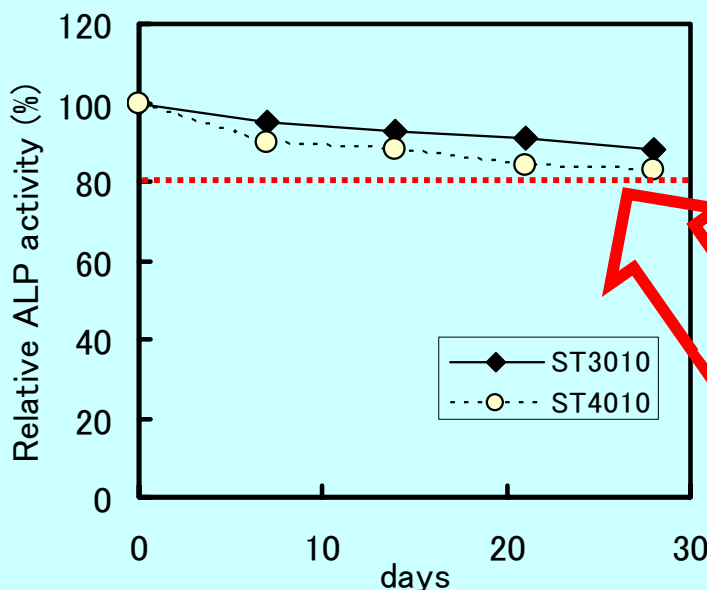

(Method)

ALP conjugate was diluted 100 times by ST3010 or ST4010.

The solutions were stored at 37 degree C for 4 weeks. The ALP activities were estimated by ALP assay kit.

0 days activity was indicated as 100%

Result shows ALP activity retained at least 80% after 4 weeks at 37°C

Description	Catalog #	volume	Price
ALP stabilizer-ST3010	51005015	100ml	¥20000 / \$191 / €154
ALP stabilizer-ST4010	51005016	100ml	¥20000 / \$191 / €154

Note: Store at 4 °C 0.09% sodium azide

 **NOF CORPORATION**

Distributor



**COSMO BIO Co., LTD.**  
Inspiration for Life Science

TOYO 2CHOME, KOTO-KU, TOKYO, 135-0016, JAPAN

http://www.cosmobio.co.jp

Phone : +81-3-5632-9617

e-mail : export@cosmobio.co.jp

FAX : +81-3-5632-9618