Code No.KAL-KR090


For research use only

Anti Phototropin 2 Polyclonal Antibody

Package Size	25 μ g (100 μ L / vial)
Format	Rabbit polyclonal antibody 0.25mg/mL
Buffer	PBS [containing 2% Block Ace as a stabilizer, 0.1% Proclin as a bacteriostat]
Storage	Store below -20°C Once thawed, store at 4°C. Repeated freeze-thaw cycles should be avoided.
Purification method	This antibody was purified from rabbit serum immunized with partial peptide of Arabidopsis PHOT2 by peptide affinity chromatography.
Working dilution	For western blotting ;0.1~0.3 μ g/mL


Western blotting

A)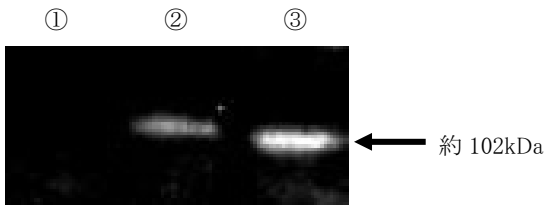

- A) Sample : Arabidopsis leaves
 ① Mutant type lacking PHOT2
 ② Hetero type
 ③ Wild type

B)

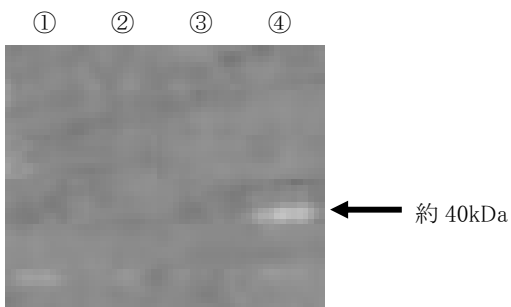- B) Sample : GST fusion protein , expressed in Escherichia coli
 ① GST-PHOT1 N terminal (without IPTG induction)
 ② GST-PHOT1 N terminal (with IPTG induction)
 ③ GST-PHOT2 N terminal (without IPTG induction)
 ④ GST-PHOT2 N terminal (with IPTG induction)

Preparation of antibodies and instruction :

Higashi T, Kagawa T, Kimura M
 Life Science and Bioengineering,
 Graduate School of Life and Environmental Science,
 University of Tsukuba

Distributor

COSMO BIO CO., LTD.
Inspiration for Life Science

TOYO 2CHOME, KOTO-KU, TOKYO, 135-0016, JAPAN

<http://www.cosmobio.co.jp>

e-mail : export@cosmobio.co.jp

Phone : +81-3-5632-9617

FAX : +81-3-5632-9618

Anti Phototropin 2 Polyclonal Antibody

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