


<u>User guide</u>




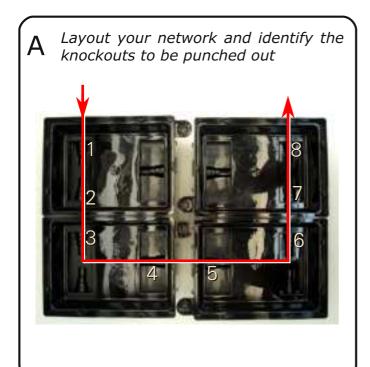

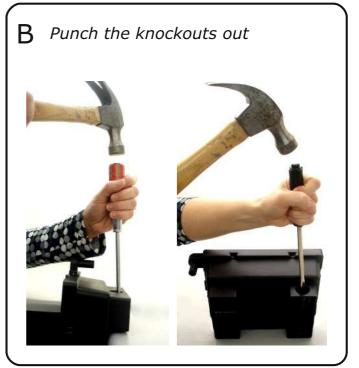

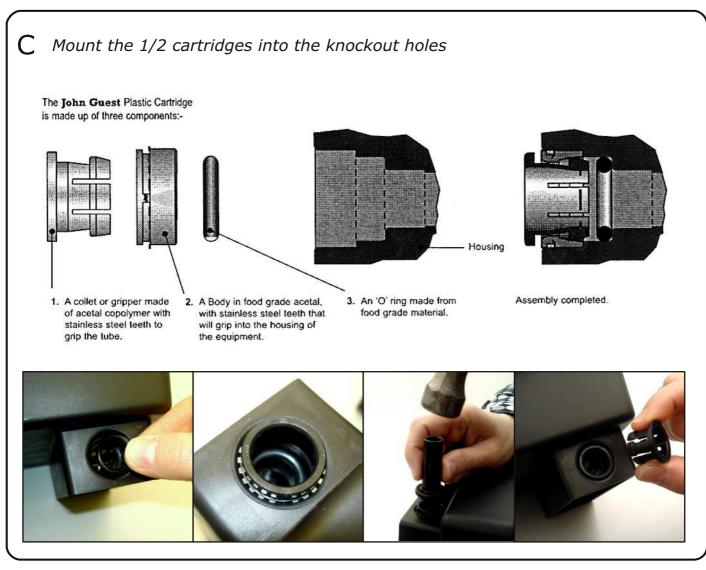

D Cut the tubing at the appropriate length and fit it into the cartridge

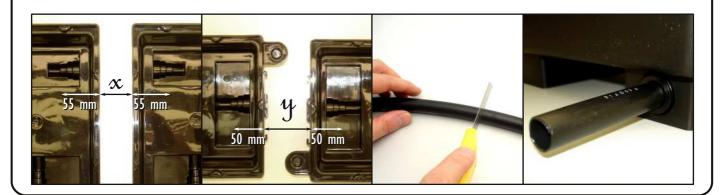

E Prepare nutrient solution from stocks




If you use FloraSeries nutrient stocks provided by Araponics, a 2000x dilution of each solution gives a good starting concentration. Nutrient replenishment and solution replacement have to be adapted to your growing conditions.

Refer to

Tocquin et al. (2003) BMC Plant Biology 2003, 3:2 http://www.biomedcentral.com/1471-2229/3/2 and references therein for more details and other formulations.

F Prepare 0.65% agar for filling the seedholders

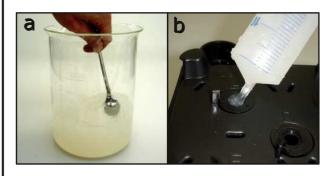

- 1. Prepare a 0.65% agar gel with double distilled water.
- 2. When solid, make it mushy (a).
- 3. Fill the seedholders when fitted (b).
- 4. Sow the seeds on the top of agar.

We wish you successful hydroponics with



