



人と科学のステキな未来へ

コスモ・バイオ株式会社

HISTOPRIME®

CatNo E6001

HistoPerm

Permanent Mounting Medium for Immunohistochemistry

Lot/Charge: See label

Storage: 24°C

Expiration Date: See label

Background

HistoPerm is a universal permanent mounting medium of tissue specimens which have been stained with histochemical and immunocytochemical methods. **HistoPerm** is a non-aqueous mounting medium offering fast drying, excellent color retention and refractivity. DAB, HistoGreen, TMB, Vector® VIP, Vector® SG, Vector® Red, Vector® NovaRED, Vector® Black und BCIP/NBT can mounted with **HistoPerm**. It is compatible with many clearing agents, including xylene and toluene.

Refraction index

1.495

Amount

100 ml **HistoPerm**

Application

HistoPerm is ready to use.

To mount tissues or cells on a slide, dispense 1 or 2 drops of **HistoPerm** onto the section. Then coverslip and allow **HistoPerm** to disperse over the entire section.

Storage

Room temperature

Note

HistoPerm contains Xylene and is flammable. Harmful by inhalation and in contact with skin. Irritating to skin.

E6001 0706271/1

AbCys S.A. - 5, rue Pierre Chausson 75010 PARIS

Tél. : 01 40 03 89 14 – Fax : 01 44 52 92 69

e-mail : infos@abcysonline.com - www.abcysonline.com