**For research use only. Not for clinical diagnosis.**

# Anti 4R-tau

## 【 Catalog No. TIP-4RT-P01】

July 1, 2024

### I. Product Overview

Tau is a microtubule-associated protein.

In adult human brain, 6 tau isoforms are expressed, which are divided into two groups, 3-repeat (3R) and 4-repeat (4R) tau isoforms. Tau assembles into insoluble, filamentous, and hyperphosphorylated inclusions in many neurodegenerative diseases, and different tau isoforms are accumulated in the diseased brains, namely, 6 tau isoforms in AD, 3R tau isoforms in Pick's disease, and 4R tau isoforms in PSP and CBD.

Therefore, isoform-specific tau antibodies are useful tools for immunohistochemical and biochemical studies of tau species in diseased brains. Tau deposited in AD is extensively deamidated at N279. This anti-4R specifically recognizes 4R tau isoforms regardless of deamidation and strongly stained tau in AD brain.

### II. Product information

<b>Product type</b>	Primary antibody
<b>Immunogen</b>	synthetic N279D 4R-tau peptide (VQIIDKKLDLSNVQSKC)
<b>Raised in</b>	Rabbit
<b>Myeloma</b>	-
<b>Clone number</b>	-
<b>Isotype</b>	-
<b>Source</b>	Anti-serum
<b>Purification</b>	-
<b>Form</b>	Liquid containing 50% glycerol and 0.05% NaN <sub>3</sub>
<b>Concentration</b>	-
<b>Volume</b>	50 uL
<b>Label</b>	Unlabeled
<b>Specificity</b>	4R-tau isoforms
<b>Cross reactivity</b>	Human, Mouse, Rat
<b>Storage</b>	Store below -20°C. (below -70°C for prolonged storage). Aliquot to avoid cycles of freeze/thaw.

### III. Technical information

#### Recommended dilutions

- Western blotting: 1/2000 - 1/5000 (Figure. A).
- Immunohistochemistry: 1:3000 - 1/20000 (Figure. B) a higher dilution is recommended for immunohistochemistry.

\*Other applications have not been tested.

\*Optimal dilutions/concentrations should be determined by the end user.

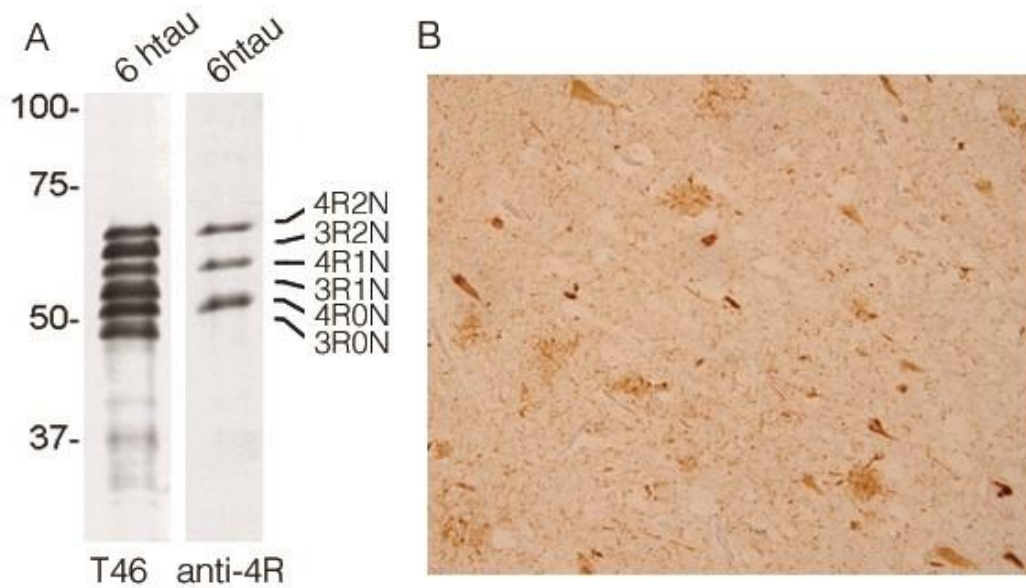
Figure A: Immuno blot analysis for 6 kinds of tau isoforms (3-repeat and 4-repeat) with anti total tau (T46) and Anti 4R (Cat No. TIP-4RT-P01)

Figure B: Immuno staining for paraffin embedded section of cerebral cortex prepared from AD brain (retrieved with proteinase K and formic acid)

#### IV. References

- 1) Hasegawa M, Watanabe S, Kondo H, et al, 3R and 4R tau isoforms in paired helical filaments in Alzheimer's disease. *Acta Neuropathol.* 2013 Nov 9
- 2) Dan A, Takahashi M, Masuda-Suzukake M, et al. Extensive deamidation at asparagine residue 279 accounts for weak immunoreactivity of tau with RD4 antibody in Alzheimer's disease brain. *Acta Neuropathol.*

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