



CEL-BSA/N^ε-(carboxyethyl) lysine-BSA

Product Description	Bovine serum albumin (BSA) (50 mg/ml) was incubated at 37°C for 24 h with pyruvate and 100 mM sodium cyanoborohydride in 0.2 M sodium phosphate buffer (pH 7.8), followed by dialysis against PBS. The CEL content (2.6 mol CEL/mol BSA) was determined by amino acid analysis.
Volume	200 ul
Concentration	1 mg/ml
Storage	Store below -20°C (below -70°C for prolonged storage). Aliquot to avoid cycles of freeze/thaw.

References

1. Nagai R., Fujiwara Y., Mera K., Yamagata K., Sakashita N., Takeya M. Immunochemical detection of N^ε-(carboxyethyl)lysine using a specific antibody. J. Immunol. Methods 332, 112-120 (2008) PMID: [18242632](#)
2. Koito W., Araki T., Horiuchi S., Nagai R. Conventional antibody against N^ε-(Carboxymethyl)lysine (CML) shows cross-reaction to N^ε-(Carboxyethyl)lysine (CEL): Immunochemical quantification of CML with a specific antibody. J Biochem. 136, 831-837 (2004) PMID: [15671494](#)
3. Nagai R., Araki T., Hayashi CM., Hayase F., Horiuchi S.. Identification of N^ε-(carboxyethyl)lysine, one of the methylglyoxal-derived AGE structures, in glucose-modified protein: mechanism for protein modification by reactive aldehydes. J Chromatogr B Analyt Technol Biomed Life Sci. 788, 75-84 (2003) PMID: [12668073](#)

Characterization

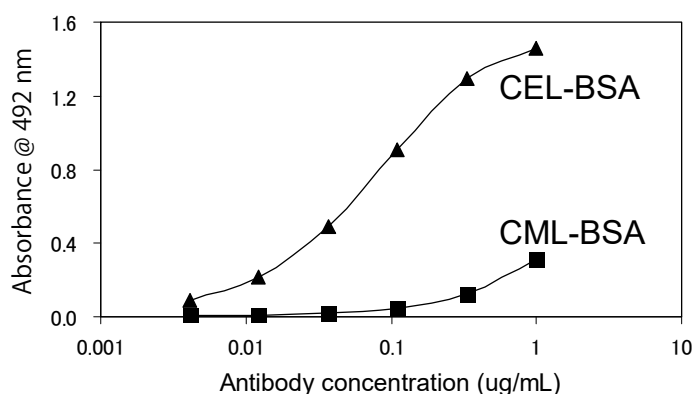


Fig.1 Immunoreactivity of CEL(CEL-SP) monoclonal antibody to CEL-BSA and CML-BSA

ELISA protocol

Coating

- 1) Distribute 100 µl / well of the sample solution (1 µg/mL in PBS) to 96 well microtiter plates (Thermo, MaxiSorp).
- 2) Incubate the plates 2h at RT or overnight at 4 degrees.
- 3) Discard the supernatant of sample solution.
- 4) Wash the plates three times with washing.buf.(PBS/0.05%Tween 20)

Blocking

- 1) Distribute 200 µl / well of 0.5% gelatin-PBS to 96 well microtiter plates
- 2) Incubate the plates 1h at RT.
- 3) Discard the the supernatant of 0.5% gelatin-PBS
- 4) Wash the plates three times with washing.buf.(PBS/0.05%Tween 20)

Primary antibody

- 1) Distribute 100 µl / well of Primary antibodies diluted with washing buf. as suggested in the APPLICATIONS to each well.
- 2) Incubate the plates 1h at RT.
- 3) Discard the supernatant of Primary antibody solution.
- 4) Wash the plates three times with washing.buf.(PBS/0.05%Tween 20)

Secondary antibody

- 1) Distribute 100 µl / well of secondary antibodies (HRP-anti mouse IgG) diluted with washing buf. as suggested in the APPLICATIONS to each well.
- 2) Incubate the plates 1h at RT.
- 3) Discard the supernatant of secondary antibody.
- 4) Wash the plates three times with washing.buf.(PBS/0.05%Tween 20)

OPD color reaction

- 1) Reaction for 2-10 minutes at RT..
- 2) Distribute 100 µL / well of 2M H₂SO₄ to each well and stop enzyme reaction.
- 3) After gentle mixing, determine the absorbance at 492 nm of each well by a spectrophotometer.

RELATED PRODUCTS:

Product Name	Quantity	Maker	Cat#
Anti N ^ε -(carboxymethyl) lysine [CML] (2G11) Monoclonal Antibody	100 ul	CAC	AGE-M01
Anti N ^ε -(carboxyethyl) lysine [CEL] (CEL-SP) Monoclonal Antibody	100 ul	CAC	AGE-M02
Anti GA-pyridine (2A2) Monoclonal Antibody	100 ul	CAC	AGE-M03
Anti N ^ω -(carboxymethyl) arginine [CMA] (3F5) Monoclonal Antibody	100 ul	CAC	AGE-M04
CML-BSA/Nε-(carboxymethyl) lysine-BSA	200 ul	CSR	AGE-GP01
CEL-BSA/Nε-(carboxyethyl) lysine-BSA	200 ul	CSR	AGE-GP02
GA-BSA/Glycolaldehyde-BSA	200 ul	CSR	AGE-GP03
Ribose-gelatin	500 ul	CSR	AGE-GP04
Mild-AGE-BSA	200 ul	CSR	AGE-GP05

For research use only, Not for diagnostic use.



COSMO BIO Co., LTD.

【JAPAN】

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU. TOKYO 135-0016, JAPAN
Phone: +81-3-5632-9610
FAX: +81-3-5632-9619
URL: <https://www.cosmobio.co.jp/>



COSMO BIO USA

【Outside Japan】

2792 Loker Ave West, Suite 101
Carlsbad, CA 92010, USA
email: info@cosmobiousa.com
Phone/FAX: (+1) 760-431-4600
URL: www.cosmobiousa.com