



ENZYME

For research use only, Not for diagnostic use.

Catalog No. UOM001**Chondroitinase ABC endo enzyme (*E.coli*)**

CAS RN 9024-13-9

EC 4.2.2.20 (Formerly EC 4.2.2.4)

Background :

Chondroitinase ABC (ChABC), an enzyme isolated and purified from *Proteus vulgaris*, a gram-negative bacillus, has the ability to degrade glycosaminoglycans such as chondroitin sulfate proteoglycans (CSPG). ChABC has been shown to degrade and inactivate CSPG in *in vivo*, and administration of the enzyme at the site of spinal cord injury accelerates recovery in many models. CSPG is known as a glycoprotein which constitutes connective tissue such as cartilage, and it shoulders the inhibitory role in nerve regeneration in the nervous system. Therefore the significance of the enzyme as a tool for studying especially in the field of neuroscience is increasing.

Product Description :

The product was purified from BL21AI *E.coli* culture, expression induced by IPTG & L-arabinose. Cells were lysed by passing through a pressure chamber in a buffer containing EDTA-free protease inhibitors. Lysate was loaded onto a 1-ml nickel-sepharose resin column in the presence of 20mM imidazole and eluted with an increasing imidazole concentration gradient in 50mM Tris-HCl + 0.5M NaCl at pH8.0. Eluted enzyme was buffer-exchanged through Sephadex G25 resin into 50mM Tris-HCl + 50mM NaCl (pH 8.0). An equal volume of glycerol was added (50% final conc.) before storing enzyme solution at -20°C.

Molecular mass (ref 1, 2): 120 ~ 145 kDa (gel filtration and sucrose gradient ultracentrifugation)

SDS-PAGE yielded 1 non-identical subunit with molecular masse of c.a. 115 kDa.

pH Optimum (ref 1): pH 8.0 (chondroitin sulfate), pH 6.8 (hyaluronic acid)**Temperature optimum (ref 1):** 37 °C**Activator (ref 2):** 0.05 M acetate**Inhibitor (ref 2):** 1 mM Zn²⁺

Specific Activity: 350mU per ml. (Activity was measured by determination of unsaturated disaccharide released from sodium chondroitin sulfate C (CSC) [Cosmo Bio Co., Ltd. CSR-NaCS-C2(ShC)3] according to the method of Morgan-Elson with N-acethyl-D-galactosamine (GalNAc) [Fuji Film Wako 019-12823] as a standard.).

Unit definition: One unit is defined as the amount of enzyme required to release of 1.0umole of unsaturated disaccharide from CSC per minute at pH 8.0, 37°C.

Other activity: essentially protease free**Precautions and Disclaimer:** This product is for Research Use Only (RUO), not for drug, household, or other uses.**Storage/Stability:** Store the product at -20 °C (DO NOT FREEZE THE ENZYME).

References:

1. Yamagata, T., et al., J. Biol. Chem., 243, 1523-1535 (1968).
2. Martinez, J.B., et al., J. Biol. Chem., 234, 2236 (1959).
3. Saito, H., et al., J. Biol. Chem., 243, 1536-1542 (1968).
4. Suzuki, S., et al., J. Biol. Chem., 243, 1543 (1968).
5. Oike, Y., et al., J. Biol. Chem., 257, 9751 (1982).

For research use only, Not for diagnostic use.



COSMO BIO Co., LTD.

【JAPAN】

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU. TOKYO 135-0016, JAPAN
Phone: +81-3-5632-9610
FAX: +81-3-5632-9619
URL: <https://www.cosmobio.co.jp/>



COSMO BIO USA

【Outside Japan】

2792 Loker Ave West, Suite 101
Carlsbad, CA 92010, USA
email: info@cosmobioussa.com
Phone/FAX: (+1) 760-431-4600
URL: www.cosmobioussa.com