## Protein Ladder

# SIMASIMA

## Instruction Manual

**COSMO BIO Co., LTD.**  
Inspiration for Life Science

### Prestained SIMASIMA-Ladderの特徴

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### Prestained SIMASIMA-Ladder/Features

- ・ The ladder is a mixture of highly purified colored recombinant proteins, and is designed for monitoring the progress of SDS-PAGE and for assessing transfer efficiency onto membranes after Western blotting.
- ・ Ready-to-use – supplied in a loading buffer for direct loading on gels; no need to boil
- ・ Use pre-stained protein ladders to determine the approximate molecular weight of proteins. For accurate molecular weight estimation of proteins, use unstained protein ladders.
- ・ Store the ladder for up to one year at 4°C.
- ・ Apparent molecular weight values for pre-stained protein markers can be different when run on different gel types. For accurate molecular weight estimation of proteins, use unstained protein ladders.

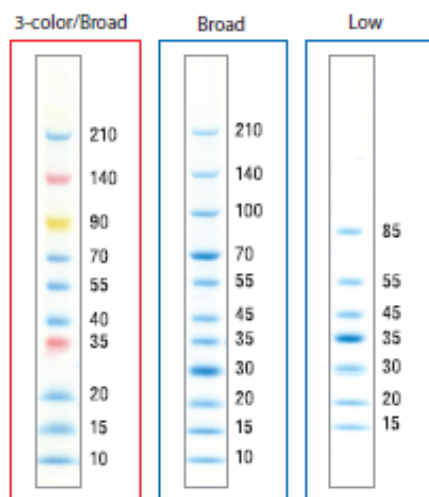
### SIMASIMA-Ladder/Features

- ・ The ladder is a mixture of highly purified recombinant 6xHis-tag fusion proteins, and is designed for precise sizing of proteins by SDS-PAGE.
- ・ Ready-to-use – supplied in a loading buffer for direct loading on gels; no need to boil
- ・ Store the ladder for up to one year at 4°C.
- ・ The molecular weight of each protein was measured by a mass spectrometry device.
- ・ Protein volume of each band when you apply 5  $\mu$ L of the ladder is 0.15  $\mu$ g (Reference band is 0.6  $\mu$ g).
- ・ Recommended load volume is 5  $\mu$ L (CBB Staining) or 0.2-0.5  $\mu$ L (Silver Staining and SYPRO Ruby Staining).

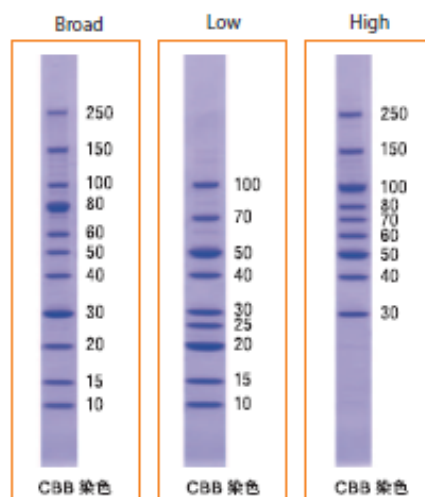
**【Note】** Prior to use, please confirm that the ladder completely dissolves.

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