# PRODUCT INFORMATION Protein A

Catalog Number: AP-01-1

Description: Pure Protein A, Alkaline Phosphatase conjugated.

Lot Number:

Protein 1 mg purified Protein A Alkaline Phosphatase / 1 ml Buffer. Concentration:

Purification Gel filtration performed after conjugation to remove free Alkaline

Procedure: Phosphatase.

Activity: 1 mg of Protein A will bind 12-15mg Human IgG.

**Buffer:** 0.01M Phosphate - 0.15M NaCl, pH 7.2 - 7.4. Contains 0.05% sodium

azide as a preservative.

Chemical Used for Conjugation:

Alkaline Phosphatase.

Storage: Store liquid refrigerated at 5-8°C in aliquots. DO NOT FREEZE!

(20-50% Glycerol has been added to prevent freezing)

Stability: The liquid material is stable for at least 1 year when stored refrigerated in

aliquots with 0.05% sodium azide added as a preservative.

Caution: Refer to the enclosed MSDS for information regarding lectins. The

aluminum seals have sharp edges and the vial itself may have cracks

which can cause lacerations. Use caution when opening the vial.

**Procedure for Use:** See reverse side.

107 Jord Amphlett Blvd.

Remarks: Protein A binds IgG from different species to varying degrees. For

details refer to Lindmark, R. et al., (1983) I. Immuno. Meth.: 61.

ABORATORIES, INC.

Tel: 650-342-3296 650-342-2648 Fax: Orders: 1-800-821-0044

(Outside CA only)

## PRODUCT INFORMATION Alkaline Phosphatase Enzyme Activity Assay

**Chemical Principle:** En

Orthophosphoric Monoester +  $H_2O \rightarrow Alcohol + H_3PO_4$ 

0.1 M Tris buffer, pH 8.2. **Assay Reagents:** BUFFER:

> ENZYME: Dilute with 0.1 M Tris Buffer.

> > Acceptable dilution: 5-20 µg/ml.

SUBSTRATE: 0.001 M p-nitrophenylphosphate (P-NPP).

Procedure:

1. Add 2.9 ml substrate to Reaction test tube and 2.9 ml to

Control test tube.

2. At time = 0, add  $100\mu$ l of diluted ENZYME to Reaction tube

and 100 µl Tris to Control tube. Mix thoroughly.

3. Measure and record optical density at 410 nm OD(410) every 15 seconds for 3 minutes, or take end point reading after 3 minutes by stopping reaction with 100µl of 5.0 M NaOH.

4. Use the OD(410) measurement to determine the rate of change

in absorbance per minute.

**Enzyme Activity** Calculations:

One unit of activity is the amount of enzyme to decompose 1 umole of P-NPP/minute at 25°C. 1.62 X 10<sup>4</sup> cm<sup>-1</sup> is the molar

absorbance of P-NPP.

OD(410) / min

OD(410) / 3min - OD(410) Control / 3 minutes

3 minutes

mg enzyme / ml reaction mixture =

[enzyme dilution]

OD(410) / min units / mg =  $1.62 \times 10^4$  ml reaction mixture

I LABORATORIES, INC.

107 North Amphlett Blvd. San Mateo, CA 94401

Tel: 650-342-3296 650-342-2648 Fax: Orders: 1-800-821-0044 (Outside CA only)

## **MATERIAL SAFETY DATA SHEET**

Effective Date: March 31, 2006 Revision 4 Page 1 of 2

#### PRODUCT IDENTIFICATION

Name: Purified proteins or biotin labeled with Horseradish Peroxidase or Alkaline

Phosphatase.

Catalog Number (s): HP-02, BA-104, BA-105, BA-108, BA-109, H-1102 to H-9000, LA-1104 to LA-9000, PA-2100 to PA-2701, AA-2100 to AA-2701, HAF-001 to HAF-2354, AAF-

001 to AAF-2354, HA-01 to HA-013, AA-01 to AA-013, HAL-1104 to HAL-

4701. AAL-1104 to AAL-4701.

Synonyms: Protein A, Avidin (egg white), Biotin, Lectins, Secondary Antibodies labeled with

Horseradish Peroxidase or Alkaline Phosphatase.

## **EMERGENCY INFORMATION**

EY Laboratories, Inc. 107 North Amphlett Blvd. San Mateo, CA 94401

**EMERGENCY PHONE:** 650 342 3296

## **HAZARDOUS COMPONENTS**

Specific protein(s) as listed on the vial label. Solutions are at a concentration generally greater than 0.5mg protein/ml. Biological activity of these labeled proteins will vary. Horseradish Peroxidase and Alkaline Phosphatase are both potent enzymes which may be harmful if ingested, inhaled, or allowed to absorb through the skin. Both enzymes are known to cause allergic responses in sensitive individuals.

## **HEALTH HAZARD INFORMATION**

EXPOSURE LIMITS: None established. The toxicological properties of these products have not been

thoroughly investigated. Care should be taken when handling any of these

**EFFECTS OF** May causes localized eye, skin, or mucous membrane irritation. Some sensitive

OVEREXPOSURE: individuals may develop a chronic allergic reaction with exposure.

Inhalation of powders and skin contact with liquids are the primary routes of **ROUTES OF EXPOSURE:** exposure. Care should be taken to avoid the formation of aerosols when

handling any of the solutions.

#### PHYSICAL CHARACTERISTICS

APPEARANCE: Powders are a light brown. Solutions will be light to dark brown. SOLUBILITY: Powders are completely soluble in many biological buffers and water.

All liquids are completely miscible in water and biological buffers.

## FIRE AND EXPLOSION HAZARDS

107 North Amphlett Blvd. San Mateo, CA 94401

**EXTINGUISHING MEDIA:** 

SPECIAL FIRE FIGHTING PRECAUTIONS:

NOTE:

Not considered to be a fire hazard.

Water spray or CO2.

None required.

Alkaline Phosphatase conjugates contain less than 0.05% sodium azide as a preservative. Azide may react with lead and copper plumbing to form explosive metal azides. Flush with copious amounts of water when disposing material in the sink.

**LABORATORIES**, INC. 650-342-3296 650-342-2648 Orders:

1-800-821-0044 (Outside CA only) MSDS for Horseradish or Alkaline Phosphatase Labeled Proteins & Biotin Continued - page 2 of 2.

REACTIVITY DATA

STABILITY: Stable. The nature of any decomposition products are not

known. They are not believed to be hazardous.

HAZARDOUS POLYMERIZATION: Will NOT occur.

INCOMPATIBILITY: None known. (Lead and copper may react with sodium

azide).

**SPILL / LEAK PROCEDURES** 

MATERIAL RELEASE / SPILL: Avoid contact with powder or liquid. Clean up spill with a

> paper towel soaked in household bleach. Do not allow solutions to dry on environmental surfaces. Wash affected area with detergent after the area has been treated with

WASTE DISPOSAL: Incinerate, autoclave, or dispose of paper waste in

> accordance with all Local, State, and Federal regulations. Due to the small quantities of material involved these products are generally not considered to be environmental hazards. All of these proteins are fully biodegradable.

### **EMERGENCY FIRST AID PROCEDURES**

May be harmful if swallowed, inhaled, or allowed to absorb through the skin. Wash contacted area with water for 15 minutes. If inhaled remove to fresh air, Report exposure to the appropriate safety official. Consult a physician if irritation occurs or if there is any indication of an allergic response, such as watering eyes, sneezing, or difficulty breathing. Any eye contact should be reported to a physician immediately

## SPECIAL HANDLING PRECAUTIONS

VENTILATION: No special ventilation is required but it is recommended to

handle these reagents in a fume hood when possible.

EYE PROTECTION: Required. Goggles or safety glasses with a side shield are

recommended.

RESPIRATORY PROTECTION: Recommended as a safety precaution, specifically when

working with powders. An approved respirator may be required for those individuals already known to be sensitive

to these materials.

PROTECTIVE GLOVES: Required when handling any of these materials.

## **SPECIAL PRECAUTIONS**

This material is for research and experimental application only. It is not intended for food, drug, household, agricultural, or cosmetic use. All materials should be handled only by technically qualified individuals experienced with working with potentially hazardous chemicals. The above information is correct to the best of our knowledge. The user should make independent decisions regarding completeness of the information, based on all sources available. EY Laboratories, Inc. shall not be held liable for any damage resulting from handling or contact with the above product.

 ${f E}\,{f Y}$  laboratories, inc.

107 North Amphlett Blvd. San Mateo, CA 94401

Tel: 650-342-3296 Fax: 650-342-2648 Orders: 1-800-821-0044 (Outside CA only)