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<http://www.fzme.com>

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TEL:045-845-2699 FAX:045-845-2889

Biological Indicator (Chlorine Dioxide)

Instruction of Use

Strain

Bacillus atrophaeus NBRC13721 (ATCC 9372)

Culture Medium

Medium: Soybean Casein Digest broth

Indicator: Phenol red

Storage

- Store under the conditions of 2-25°C, 20-80% RH, away from light. (Store the culture medium under the conditions of 2-10°C)
- Do not store these indicators near sterilants or other chemicals.

Before using

- Please check expiration date before using.
- Please check the integrity before using.
- Let the indicators return to room temperature before using, in the case of refrigerated storage.

Disposal

- Dispose of any positive cultures or expired indicators by autoclaving at 121°C for not less than 30 minutes.

Protocol

- 1 Place spore strip bags in the most challenging areas.
- 2 Following exposure to CD gas, remove the bags as soon as the area has been aerated. Using strict aseptic clean bench, transfer each strip to a tube of culture medium.
- 3 (Use ACE mini Incubator Model H8200) Incubate the strips at the temperature of 30-37°C for 48 hours or 7 days.
- 4 As a positive growth control, transfer a non-sterilized strip to culture medium and incubate meanwhile. Positive control should be included in each test.

Interpretation

Negative ⇒ No color change ⇒ Adequate sterilation

Positive ⇒ A color change toward yellow ⇒ Bacterial growth

COSMO BIO CO., LTD.

TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
KOTO-KU, TOKYO 135-0016, JAPAN
TEL : (81)3-5632-9617
FAX : (81)3-5632-9618
e-mail : export@cosmobio.co.jp
URL : www.cosmobio.com

www.cosmobio.com