Product data sheet

# COSMEDIUM 005

## Serum-free liquid medium for hybridoma cell

Cat. No. COS-005

Published: 09/12/2022

### Applications

This product is a serum-free liquid medium for mouse and human hybridoma cultures that has been developed for application in mass-production of antibodies.

### Features

This product enables high-density culture of hybridomas without depletion of nutrients, even if the medium is not replaced for about one week.

### Specification

- Packaging: 1000mL
- Storage conditions: 4°C (in a cool and dark place)

Inspection Items	Specifications
Appearance	Yellow-orange liquid, no precipitation
Sterility* <sup>1</sup>	Negative
Mycoplasma* <sup>2</sup>	Negative
pH* <sup>3</sup>	7.10±0.20
Endotoxin* <sup>4</sup>	0.3 EU/mL or less
Osmolarity (mOsm/kg)* <sup>5</sup>	300±20 mOsm/kg
Cell proliferative performance* <sup>6</sup>	More than 75% of control
Cell proliferative performance* <sup>7</sup>	More than 75% of control

\*1 Culture on thioglycolic acid medium I and SCD medium for sterility test (after 14 days).

\*2 Detection by polymerase chain reaction (PCR).

\*3 Measured after incubation at 37°C for at least 12 hours under 5% CO<sub>2</sub> gas.

\*4 Measurement by kinetic colorimetric method.

\*5 Measurement by freezing point depression method.

\*6 Proliferation performance against hybridoma cells for the previous lot.

\*7 Proliferation performance against lymphocyte cells for the previous lot.

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