[Cell Freezing Medium] (Chemically defined)
[Cell Freezing Medium]

COS banker COS banker II

Catalog No. KOJ-COS-CFM01

Catalog No. KOJ-COS-CFM02

COS banker is a **one-step culture cell storage solution**. Suspend culture cells in COS banker and store at -80°C for long term storage (it is not necessary to store in liquid nitrogen). This is a serum-free solution with no animal products added.

Features

- ▶ Ready-to-use
- ▶ No dilutions required
- ▶ Store in -80°C freezer
- ▶ Rapid and long term storage at -80°C
- ▶ No serum
- ▶ No animal products

Instructions




1. Collect culture cells (during logarithmic growth stage) and place 5×10^5 - 1×10^7 cells in a cryovial. Centrifuge for 5 minutes at 1,200 rpm.
2. Add 1 ml of COS banker to cell pellet and vigorously mix, then freeze and store at -80°C.
3. If liquid Nitrogen storage is desired, freeze cells at -80°C first and then transfer to liquid Nitrogen.
4. To thaw, defrost in 37°C water bath quickly.

Note: It is ideal to use a logarithmic growth stage culture in order to increase the survival rate when thawed.

Contents and Usage

Product Name	COS banker	COS banker II
Product Code	KOJ- COS-CFM01	KOJ- COS-CFM02
Size	120 mL	120 mL
Specification	Contains DMSO No serum	Contains DMSO Contains Human Serum Albumin No serum
Storage	Store at 2-8°C of dark place	Store at 2-8°C of dark place
Expiry Date	2 years from date of manufacture	2 years from date of manufacture
Application Tested Cell	Daudi, HeLa S3, K562, Raji, MOLT-4, CTLL-2, NIH/3T3, PC-12, CHO-K1	Activated lymphocytes derived from human peripheral blood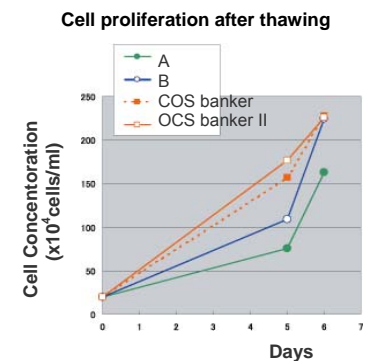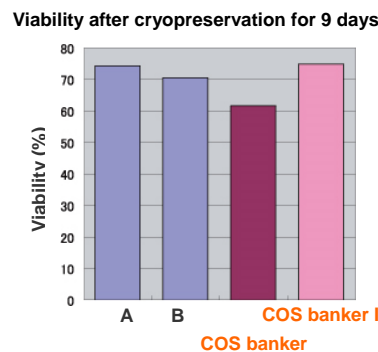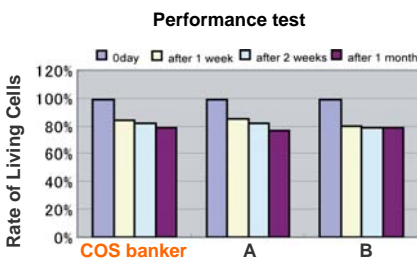

Quality Controls

Endotoxin-free tested / Bacteria free tested

Warnings

- ▶ Research use only
- ▶ Not for diagnostic and/or clinical use

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[Cell Freezing Medium] (Chemical defined)

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COSMO BIO CO., LTD.