# Gelatin-Zymography Precast GEL (Cosmo Bio type)

Catalog No. AK48, AK50

April 2, 2020

## Introduction

Matrix metalloproteinases (MMPs) belong to the family of metalloproteinases, which consists of at least 20 members, and known to be involved in the metabolism of extracellular matrix proteins. MMPs are widely detected by zymography. The Gelatin-Zymography Kit (Cat.No.AK47) provides an easy system of the electrophoresis for zymography. This product is used for detecting ProMMP-2, MMP-2 and ProMMP-9 and MMP-9 in blood, body fluid, secretion, cell lysate, cell culture medium, and other samples.

## Components/Storage

Catalog No. / Component	Quantity	Storage
AK48 / Precast gel for 12 samples	5 pieces	4°C
AK50 / Precast gel for 12 samples	10 pieces	4°C

The size of the glass plate gel is 100(W)×100(H) and the thickness is 1mm.

## References

- (1) Miyazaki, K., Ohta, Y., Nagai, M., Morimoto, N., Kurata, T., Takehisa, Y., Ikeda, Y., Matsuura, T., Abe, K. Disruption of Neurovascular Unit Prior to Motor Neuron Degeneration in Amyotrophic Lateral Sclerosis. *J. Neurosci. Res.* 89, 718-728 (2011)
- (2) Fujiwara, M., Kashima, T. G., Kunita, A., Kii, I., Komura, D., Grigoriadis, A. E., Kudo, A., Aburatani, H., Fukayama, M. Stable Knockdown of S100A4 Suppresses Cell Migration and Metastasis of Osteosarcoma. *Tumour Biol.*
- (3) Aoki, T., Kataoka, H., Ishibashi, R., Nozaki, K., Hashimoto, N. Simvastatin Suppresses the Progression of Experimentally Induced Cerebral Aneurysms in Rats. *Stroke.* 39, 1276-85 (2008)
- (4) Kuramochi, D., Unoki, H., Bujo, H., Kubota, Y., Jiang, M., Rikihisa, N., Udagawa, A., Yoshimoto, S., Ichinose, M., Saito, Y. Matrix Metalloproteinase 2 Improves The Transplanted Adipocyte Survival in Mice. *Eur. J. Clin. Invest.* 38, 752-759 (2008)
- (5) Fujimoto, M., Takagi, Y., Aoki, T., Hayase, M., Marumo, T., Gomi, M., Nishimura, M., Kataoka, H., Hashimoto, N., Nozaki, K. Tissue Inhibitor of Metalloproteinases Protect Blood-brain Barrier Disruption in Focal Cerebral Ischemia. *J. Cereb. Blood Flow Metab.* 28, 1674-1685 (2008)

## **Notice to Purchaser**

This product is to be used for Research Purposes Only. It is not to be used for Drug or Diagnostic Purposes, nor is it intended for Human Use. Cosmo Bio products may not be resold, modified for resale, or used to manufacture commercial products without the express written consent of Cosmo Bio Co., Ltd.

Except as otherwise expressly set forth in this use manual, Cosmo Bio does not make any representation or warranties or conditions of any kind, either express or implied, with respect to the products, or information disclosed hereunder, including, but not limited to, the implied warranties of merchantability, fit for a particular purpose, or noninfringement of the intellectual property rights of third parties.

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COSMO BIO Co., LTD.