Osteoclast Culture Medium (for human)

Catalog No. OSCMHB (Japan)
Catalog No. PMC-OSCMHB (Outside Japan)

Edited Dec 24, 2025

Please read this manual carefully before use. This product is for research use only.

Product Overview

Bone metabolism is regulated by a balance between bone formation (osteogenesis) and bone resorption. Previous studies have demonstrated that bone marrow cells can differentiate into osteoclasts in the presence of M-CSF (macrophage colony-stimulating factor) and RANKL (receptor activator of NF- κ B ligand).

Product Information

This product is a culture medium containing lot-tested serum, RANKL (receptor activator of NF- κ B ligand), M-CSF (macrophage colony-stimulating factor), and antibiotics.

This medium is designed to support the differentiation of human bone marrow cells into osteoclasts. Osteoclast cells (human) can be purchased separately (Cat. No. OSC15C (Japan) Cat. No. PMC-OSC15C (Outside Japan)).

Storage Conditions

This product is shipped in a frozen state. Upon receipt, store immediately at $-20\text{ }^{\circ}\text{C}$ or below. The expiration date indicated on the product label applies to the unopened product stored at $-20\text{ }^{\circ}\text{C}$ or below. After thawing, store at $4\text{ }^{\circ}\text{C}$ and use as soon as possible. The product may be refrozen once only. After thawing, it is recommended to aliquot into appropriate volumes for several uses and store frozen. Repeated freeze-thaw cycles may significantly compromise product quality.

Instructions for Use

Remove the product from the freezer and allow it to thaw slowly in a refrigerator. Do not warm the product prior to use. When required ready to use, aliquot the necessary volume and allow it to warm to room temperature before using. Prolonged storage at room temperature or repeated warming to room temperature may significantly compromise product quality. Use this product under aseptic conditions.

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