

## Instructions For Use ACA-IFU

Rev. Date: 10/31/03

Revision: 3

Page 1 of 1

P.O. Box 3286 - Logan, Utah 84323, U.S.A. - Tel. (800) 729-8350 - Fax (435) 755-0015 - [www.scytek.com](http://www.scytek.com)

# Peroxide Block

**Description:** Peroxide Block has been developed to use with immunolabeling techniques for the reduction of nonspecific background staining do to endogenous peroxidase.

**Appearance:** Clear liquid

**Uses:** Immunohistochemical procedures.

**Limitations:** Do not use past expiration date.

<b>Availability:</b>	<u>REF</u>	<u>Volume</u>
	ACA125	125 ml
	ACA500	500 ml
	ACA999	1000 ml

**Storage:** Store Peroxide Block at 2-8°C. Product is stable for 18 months from date of manufacture.

### Procedure:

1. Incubate tissue section for 10 minutes at either room temperature or 37° C prior to application of the primary antibody. After incubation, rinse once in buffer.
2. \*\*\* For bulk staining, pour Peroxide Block in a covered staining tray and dip slides for 5 minutes. Replace with fresh reagent after 5-10 uses.

**Precautions:** Do not pipette reagent by mouth.  
Avoid contact with skin and eyes.  
Wash after use.  
Observe all federal, state and local environmental regulations regarding disposal.

Storage: 2° C



8° C



ScyTek Laboratories, Inc.  
205 South 600 West  
Logan, UT 84321  
435-755-9848  
U.S.A.



Authorized Representative in Europe



(Regulatory affairs only)  
EmergoEurope (31)(0) 70 345-8570  
Molsstraat 15  
2513 BH Hague, The Netherlands