

Instructions For Use HMM-IFU

Rev. Date: 10/28/03

Revision: 2

Page 1 of 1

P.O. Box 3286 - Logan, Utah 84323, U.S.A. - Tel. (800) 729-8350 - Fax (435) 755-0015 - www.scytek.com

Mayer's Hematoxylin (Lillie's Modification)

Description: Mayer's Hematoxylin (Lillie's Modification) is a stain with histological applications.

Results: Nuclei are blue.

Availability:

| <u>Item #</u> | <u>Volume</u> |
|---------------|---------------|
| HMM500 | 500 ml |
| HMM999 | 1 Liter |
| HMM010 | 10 Liters |

Procedure:

1. Re-hydrate tissue as usual.
2. Stain nuclei 5 minutes with hematoxylin.
3. Place in Bluing Reagent (BRT500) until stain is blue.
4. Rinse in DI water.
5. Dehydrate, clear, and mount in permanent mounting medium.